


Water drainage for / Wasserablauf für 889 IC Sample Center

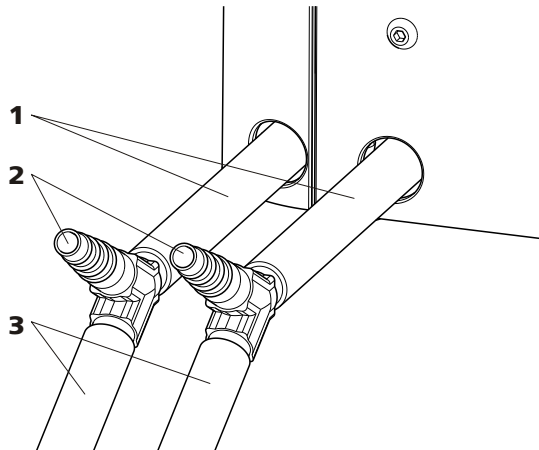

- 1 Drainage tubing to the Y connector / Ablaufschlauch zum Y-Verbinder (6.1816.080)**
- 2 Y connector for tubing 6-9 mm ID / Y-Verbinder für Schlauch ID 6-9 mm (6.1807.010)**
- 3 Drainage tubing to the waste container / Ablaufschlauch zum Abfallbehälter (6.1816.080)**

EN

Mounting the water drainage

To prevent liquid from backing up in the the water drainage tubing a Y connector must be attached for aeration.

1. Cut the drainage tubings [1] to a length of approx. 10 cm and attach to each drain nozzle.
2. Attach the Y connectors [2] to the drainage tubings.
3. Attach the drainage tubings [3] to the lower outlets of the Y connectors [3]. The length of the tubings should be selected in accordance with the height of the waste container.



Note

The waste container must be at a lower level than the drain nozzles of the 889 IC Sample Center.

DE

Wasserablauf montieren

Damit sich im Wasserablauf-Schlauch kein Rückstau bildet, ist ein Y-Verbinder zur Belüftung anzubringen.

1. Die Ablaufschläuche [1] auf ca. 10 cm zuschneiden und jeweils auf die Ablaufstutzen aufstecken.
2. Die Y-Verbinder [2] auf die Ablaufschläuche aufstecken.
3. Die Ablaufschläuche [3] auf die unteren Abgänge der Y-Verbinder [3] aufstecken. Die Länge der Schläuche richtet sich entsprechend der Höhe zum Abfallbehälter.



Hinweis

Das Niveau des Abfallbehälters muss tiefer als die Ablaufstutzen des 889 IC Sample Center sein.