850 Professional IC



AnCat – non-suppressed – 2.850.3000

Manual 8.850.8054EN / 2019-11-29





Metrohm AG CH-9100 Herisau Switzerland Phone +41 71 353 85 85 Fax +41 71 353 89 01 info@metrohm.com www.metrohm.com

850 Professional IC

AnCat – non-suppressed – 2.850.3000

Manual

Technical Communication Metrohm AG CH-9100 Herisau techcom@metrohm.com

This documentation is protected by copyright. All rights reserved.

This documentation has been prepared with great care. However, errors can never be entirely ruled out. Please send comments regarding possible errors to the address above.

Table of contents

Table of contents

1	Introdu	ctio	n	1
	1	.1	Instrument description	. 1
	1	.2	Intended use	. 1
		.3 .3.1	About the documentation Symbols and conventions	
	1 1 1 1	.4.1 .4.2 .4.3 .4.4 .4.5	Safety instructions General notes on safety Electrical safety Tubing and capillary connections Flammable solvents and chemicals Recycling and disposal	3 3 . 4 4
2	Overvie	ew o	f the instrument	6
	2	.1	Front	. 6
	2	.2	Rear	. 8
3	Installa	tion		10
	3	.1	About this chapter	10
	3	.2	Initial installation	10
	3	.3	Installation diagram	12
	3	. 4 .4.1 .4.2 .4.3	Setting up the instrument Packaging Checks Location	14 14
	3	.5	Capillary connections in the IC system	15
	3 3 3 3	.6.1 .6.2 .6.3 .6.4 .6.5	Rear of the instrument Rollers and handle Positioning and connecting the detector Transport locking screws Leak sensor Drainage tubings	17 20 20 20
	3	.7	Capillary and cable feed-throughs	23
		.8 .8.1	Eluent	
	3	.9	Eluent degasser	30
	3	. 10 .10.1 .10.2	High pressure pump Capillary connections high pressure pump/purge valve Deaerating the high pressure pump	32

		3.11	Inline filter	. 36
		3.12	Pulsation absorber	37
		3.13	Sample degasser	. 38
		3.14 3.14.1 3.14.2 3.14.3	Injection valve Connecting the injection valve Mode of operation of the injection valve Selecting the sample loop	. 40 . 41
		3.15	Column thermostat	42
		3.16	Conductivity detector	. 45
		3.17	Connecting the instrument to a computer	. 47
		3.18	Connecting the instrument to the power grid	
		3.19	Guard column	49
		3.20	Separation column	50
4	Start-	up		53
		4.1	Initial start-up	. 53
		4.2	Conditioning	. 54
5	Opera	ition a	and maintenance	56
		5.1 5.1.1 5.1.2 5.1.3 5.1.4	General notes Care	. 56 56 57
		5.2 5.2.1	Capillary connections Operation	
		5.3	Door	. 58
		5.4 5.4.1 5.4.2	Eluent Production Operation	58
		5.5 5.5.1 5.5.2	High pressure pump Protection Maintenance	59
		5.6 5.6.1	Inline filter	
		5.7	Inline sample preparation	. 72
		5.8	Rinsing the sample path	72
		5.9 5.9.1	Sample degasser Operation	

Table of contents

	5.10 5.10.1	Injection valve	
	5.11	Conductivity detector	
	5.11.1		
	5.12 5.12.1 5.12.2 5.12.3 5.12.4	Protection	75 75 76
6	Troublesho	ooting	77
	6.1	Problems and their solutions	77
7	Technical s	pecifications	81
	7.1	Reference conditions	81
	7.2	Device	81
	7.3	Leak sensor	81
	7.4	Ambient conditions	81
	7.5	Housing	82
	7.6	Eluent degasser	82
	7.7	High pressure pump	82
	7.8	Sample degasser	83
	7.9	Injection valve	83
	7.10	Column thermostat	84
	7.11	Conductivity measuring system	84
	7.12	Power connection	85
	7.13	Interfaces	85
	7.14	Weight	86
8	Accessorie	s	87
	Index		88

Table of figures

Table of figures

Figure 1	Front 850 Professional IC – AnCat – non-suppressed	. 6
Figure 2	Rear 850 Professional IC AnCat – non-suppressed	. 8
Figure 3	Installation diagram	13
Figure 4	Connection of capillaries with pressure screws	15
Figure 5	Rollers and handle	18
Figure 6	Handle as MPak holder	
Figure 7	Connection of the leak sensor on the rear of the instrument	21
Figure 8	Drainage tubings	
Figure 9	Capillary feed-throughs on the door	
Figure 10	Capillary feed-throughs base tray/flask holder	25
Figure 11	Installing the eluent bottle cap	27
Figure 12	Mounting the aspiration filter	
Figure 13	Installing the tubing weighting and aspiration filter	28
Figure 14	Eluent aspiration tubing fully equipped	28
Figure 15	Eluent bottle – connected	29
Figure 16	Eluent degasser	
Figure 17	Capillary connections high pressure pump/purge valve	32
Figure 18	High pressure pump – Connect inlet	33
Figure 19	Deaerate the high pressure pump	
Figure 20	Connecting the inline filter	
Figure 21	Pulsation absorber – Connection	38
Figure 22	Sample degasser	39
Figure 23	Injection valve – connected	
Figure 24	Injection valve – Positions	41
Figure 25	Column thermostat	43
Figure 26	Conductivity detector front	45
Figure 27	Conductivity detector rear	
Figure 28	Connection detector – separation column	47
Figure 29	Pump head – removing the piston	61
Figure 30	Components of the piston cartridge	62
Figure 31	Tool for piston seal	63
Figure 32	Removing the piston seal	64
Figure 33	Inserting the piston seal into the tool	64
Figure 34	Inserting the piston seal into the pump head	65
Figure 35	Removing valves	66
Figure 36	Dismantling valve	
Figure 37	Components of the inlet valve and outlet valve	68
Figure 38	Change filters (of the inline filter)	70

1 Introduction

1 Introduction

1.1 Instrument description

The instrument **850 Professional IC – AnCat – non-suppressed** (2.850.3000) is one of the model versions of the Professional IC line of instruments manufactured by the Metrohm Company. The Professional IC line of instruments is distinguished by

- the intelligence of its components, which are able to monitor and optimize all functions and to provide documentation according to FDA requirements.
- its compactness.
- its flexibility. A suitable model version exists for every application.
 Individual instruments can be converted, expanded or modified to create a different instrument as needed.
- its **transparency**. All components are easily accessible and arranged in a clear manner.
- its **safety**. Chemicals and electronics are separated and a leak sensor is integrated in the wet end.
- its environmental compatibility.
- its low noise emission.

The instrument is operated with **MagIC Net** software. It is connected via a USB connection to a PC on which MagIC Net is installed. The software automatically recognizes the instrument and checks its functional readiness. MagIC Net controls and monitors the instrument, evaluates the measured data and administers it in a database. The operation of MagIC Net is described in the online help or in the tutorial for MagIC Net.

1.2 Intended use

The instrument **850 Professional IC – AnCat – non-suppressed** is used for the ion chromatographic determination of cations and anions without suppression. The two channels can be operated in parallel or completely independently of one another. Thanks to the two high pressure pumps, the instrument is also suitable for applications with high pressure gradients.

The present instrument is suitable for processing chemicals and flammable samples. The usage of the 850 Professional IC – AnCat – non-suppressed therefore requires that the user has basic knowledge and experience in the handling of toxic and caustic substances. Knowledge with respect to

1.3 About the documentation

the application of the fire prevention measures prescribed for laboratories is also mandatory.

1.3 About the documentation



CAUTION

Please read through this documentation carefully before putting the instrument into operation. The documentation contains information and warnings which have to be followed by the user in order to ensure safe operation of the instrument.

1.3.1 Symbols and conventions

The following symbols and styles are used in this documentation:

(5- 12)	Cross-reference to figure legend		
	The first number refers to the figure number, the second to the instrument part in the figure.		
1	Instruction step		
	Carry out these steps in the sequence shown.		
	Warning		
/ :	This symbol draws attention to a possible life hazard or risk of injury.		
	Warning		
7	This symbol draws attention to a possible hazard due to electrical current.		
	Warning		
<u></u>	This symbol draws attention to a possible hazard due to heat or hot instrument parts.		
	Warning		
	This symbol draws attention to a possible biological hazard.		
	Caution		
	This symbol draws attention to a possible damage of instruments or instrument parts.		
•	Note		
	This symbol marks additional information and tips.		

1 Introduction

1.4 Safety instructions

1.4.1 General notes on safety



WARNING

Operate this instrument only according to the information contained in this documentation.

This instrument left the factory in a flawless state in terms of technical safety. To maintain this state and ensure non-hazardous operation of the instrument, the following instructions must be observed carefully.

1.4.2 Electrical safety

The electrical safety when working with the instrument is ensured as part of the international standard IEC 61010.



WARNING

Only personnel qualified by Metrohm are authorized to carry out service work on electronic components.



WARNING

Never open the housing of the instrument. The instrument could be damaged by this. There is also a risk of serious injury if live components are touched.

There are no parts inside the housing which can be serviced or replaced by the user.

Supply voltage



WARNING

An incorrect supply voltage can damage the instrument.

Only operate this instrument with a supply voltage specified for it (see rear panel of the instrument).

1.4 Safety instructions

Protection against electrostatic charges



WARNING

Electronic components are sensitive to electrostatic charges and can be destroyed by discharges.

Do not fail to pull the power cord out of the power socket before you set up or disconnect electrical plug connections at the rear of the instrument.

1.4.3 Tubing and capillary connections



CAUTION

Leaks in tubing and capillary connections are a safety risk. Tighten all connections well by hand. Avoid applying excessive force to tubing connections. Damaged tubing ends lead to leakage. Appropriate tools can be used to loosen connections.

Check the connections regularly for leakage. If the instrument is used mainly in unattended operation, then weekly inspections are mandatory.

1.4.4 Flammable solvents and chemicals



WARNING

All relevant safety measures are to be observed when working with flammable solvents and chemicals.

- Set up the instrument in a well-ventilated location (e.g. fume cupboard).
- Keep all sources of flame far from the workplace.
- Clean up spilled liquids and solids immediately.
- Follow the safety instructions of the chemical manufacturer.

1.4.5 Recycling and disposal



This product is covered by European Directive 2012/19/EU, WEEE – Waste Electrical and Electronic Equipment.

The correct disposal of your old instrument will help to prevent negative effects on the environment and public health.

1 Introduction

More details about the disposal of your old instrument can be obtained from your local authorities, from waste disposal companies or from your local dealer.

2.1 Front

2 Overview of the instrument

2.1 Front

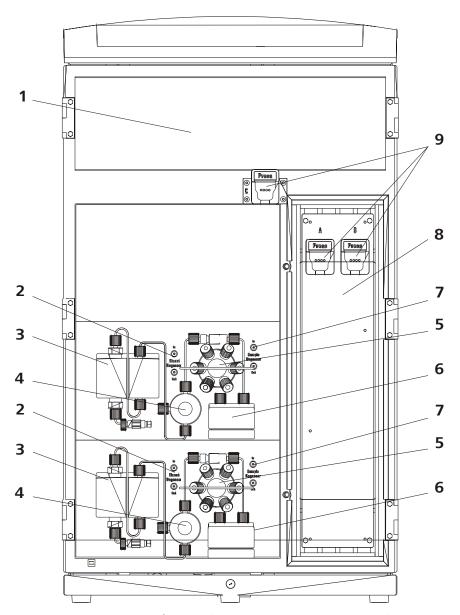


Figure 1 Front 850 Professional IC – AnCat – non-suppressed

1 Detector chamber

Space for one conductivity detector each for the two channels.

2 Eluent degasser

One eluent degasser each for the two channels.

See Chapter 3.9, Page 30

2 Overview of the instrument

3 High pressure pump

One high pressure pump each for the two channels.

See Chapter 3.10, Page 32

5 Injection valve

One injection valve each for the two channels.

See Chapter 3.14, Page 40

7 Sample degasser

One sample degasser each for the two channels. Utilization optional.

See Chapter 3.13, Page 38

9 Column holder

For two separation columns in the column thermostat and one additional separation column outside the column thermostat.

See Chapter 3.20, Page 50

4 Purge valve

One purge valve each for the two channels. See Chapter 3.10.1, page 32

6 Pulsation damper

One pulsation damper each for the two channels.

See Chapter 3.12, Page 37

8 Column thermostat

See Chapter 3.15, Page 42.

It is conventional to refer to the channel with the upper high pressure pump and the left-hand conductivity detector as the **anion channel**, and the channel with the lower high pressure pump and the right-hand conductivity detector as the **cation channel**.

2.2 Rear

2.2 Rear

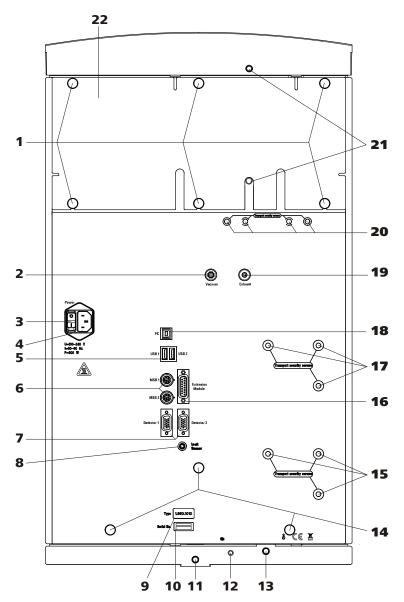


Figure 2 Rear 850 Professional IC AnCat – non-suppressed

1 Knurled screws

For fastening the rear panel (2-22) and handle (5-2).

3 Mains switch

For switching the instrument on and off.

I = ON

0 = OFF

2 Vacuum connection

For connecting further degassing chambers in extension modules.

Labeled with Vacuum.

4 Mains connection socket

For connecting the mains cable.

2 Overview of the instrument

5 USB connectors

2 USB connectors (labeled with *USB 1* and *USB 2*).

7 Detector connectors

2 detector connectors (labeled with *Detector 1* and *Detector 2* for connecting Metrohm detectors.

9 Instrument type

11 Drainage tubing connection

For connecting a drainage tubing 6.1816.020 (8-8).

13 Drainage tubing connection

For connecting a drainage tubing 6.1816.020 (8-**9**).

15 Transport locking screws

For securing the lower high pressure pump (17-4) when transporting the instrument (only necessary for instruments with two high pressure pumps).

17 Transport locking screws

For securing the high pressure pump (17-4) when transporting the instrument.

19 Waste air opening

For extracting the air from the vacuum chamber.

Labeled with Exhaust.

21 Drainage tubing connection

For connecting a drainage tubing 6.1816.020 (8-**1**).

6 MSB connectors

2 MSB connectors (labeled with MSB 1 and MSB 2) for connecting MSB devices

Caution: The 850 **must** be switched off when connecting a device.

MSB = Metrohm Serial Bus.

8 Leak sensor connection socket

For connecting the leak sensor connector plug (7-**2**).

10 Serial number

12 Leak sensor connection cable

For connecting the leak sensor.

14 Knurled screws

For fastening the rollers.

16 Extension module connection

For connecting an extension module (labeled with *Extension Module*).

18 PC connection socket

For connecting the instrument to the computer with the USB cable 6.2151.020.

20 Transport locking screws

For securing the vacuum pump when transporting the instrument.

22 Rear panel

Removable. Access to the detector chamber.



CAUTION

When connecting an instrument to the MSB connector (2-6) you must switch off the 850 Professional IC.

3.1 About this chapter

3 Installation

3.1 About this chapter

The Installation chapter contains

- this overview
- a brief set of instructions for the initial installation of the 850 Professional IC AnCat non-suppressed (see chapter 3.2, page 10). At each step you will find cross-references to more detailed installation instructions for individual components, should you require such aids.
- an installation diagram (see chapter 3.3, page 12), showing a completely installed instrument.
- several chapters with detailed installation instructions for all components, including those that are already installed at the time the instrument is delivered.

3.2 Initial installation



NOTICE

A number of the capillaries is already connected at the time the instrument is delivered.

The following work steps must still be carried out:

Installing 850 Professional IC - AnCat - non-suppressed



NOTICE

Each of the steps numbered 3, 4, and 7 are to be carried out for the anion and cation channels, respectively.

1 Setting up the instrument

See Chapter 3.4, Page 14.

- 2 Installations on the rear of the instrument
 - Remove handle and rollers (see chapter 3.6.1, page 17).

• Place the detectors in the instrument and connect them (see chapter 3.6.2, page 20).

- Remove transport locking screws (see "Removing transport locking screws", page 20).
- Connect the leak sensor (see "Connecting the leak sensor ", page 20).
- Connect drainage tubings (see "Installing drainage tubings", page 23).

3 Installing the eluent path

The following steps need to be carried out once for each of the two channels.

- Equip the eluent aspiration tubing and connect with the eluent bottle (see chapter 3.8.1, page 26).
- At the end of the pre-installed column inlet capillaries (3-2) and, respectively (3-5), connect one 6.2744.040 coupling each, using a 6.2744.014 pressure screw, instead of the columns.
- Connect detector inlet capillaries (3-3) and, respectively (3-6), to the 6.2744.014 coupling, using a 6.2744.040 pressure screw.

4 Installing the sample path



NOTICE

The sample degasser does not have to be connected. We recommend the usage of the sample degasser only if the sample matrix requires it.

The following steps need to be carried out once for each of the two channels.

- Connect the 6.1803.040 sample aspirating capillary (3-8) connected at the sample input of the injection valve to the output of the sample degasser with the aid of a 6.2744.090 long pressure screw (see chapter 3.13, page 38).
- Connect a part of the 6.1803.040 sample aspirating capillary with a 6.2744.090 long pressure screw to the input of the sample degasser. Guide the other end out of the instrument through a capillary feed-through.
- In the cation channel, replace the 6.1825.210 sample loop with the 6.1825.220 sample loop (see "", page 41).

5 Connecting the instrument

• Connect the instrument to the PC.

3.3 Installation diagram

• Connect instrument to mains supply .

6 Initial start-up

(see chapter 4.1, page 53).

- Switch on the PC and start MagIC Net.
- Switch on and prepare the instrument.
- Deaerate the high pressure pump.
- Rinse the instrument without column(s).

7 Installing the columns

The following steps need to be carried out once for each of the two channels:

- Remove the 6.2744.040 coupling between column inlet capillaries and detector inlet capillaries.
- Install the guard column (if used) (see "Connecting and rinsing the guard column", page 50).
- Install the separation column (see "Connecting and rinsing the separation column", page 51).
- Connect the detector inlet capillary (see "Connect the detector input capillary to the separation column", page 46).

8 Conditioning the instrument

(see chapter 4.2, page 54).

The instrument is now ready for measuring samples.

3.3 Installation diagram

Figure 3, page 13 shows the capillary connections of a system with a cation channel and an anion channel.

The two channels can be operated in parallel or completely independently of one another.

The arrangement of the modules in the diagram corresponds to the front view of the instrument. Liquid containers (eluent bottle, sample vessel, waste container, auxiliary solutions container) and guard column (see chapter 3.19, page 49) are not shown in the diagram.

A few of the capillaries are already pre-installed at the time the instrument is delivered. Capillaries on which nothing needs to be done at the time of initial installation are not numbered in the diagram.

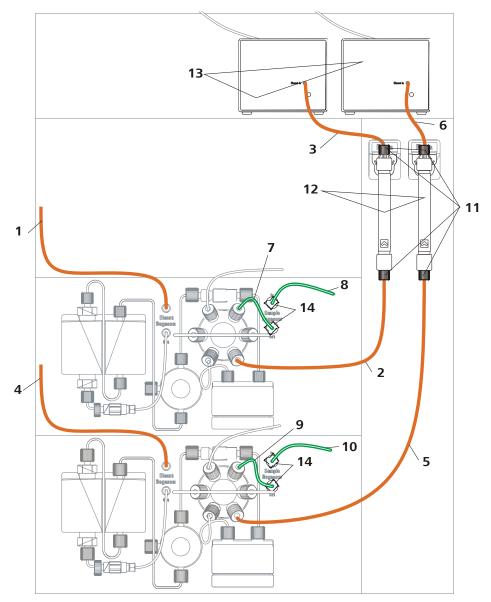


Figure 3 Installation diagram

- **1 Eluent aspiration tubing 6.1834.080** Connected to the eluent degasser.
- **3** Detector inlet capillary
- **5** Column inlet capillary

 Connected to the injection valve and threaded into the capillary recesses of the column thermostat.
- 2 Column inlet capillary
 Connected to the injection valve and
 threaded into the capillary recesses of the
 column thermostat.
- **Eluent aspiration tubing 6.1834.080**Connected to the eluent degasser.
- **6** Detector inlet capillary

7 Sample aspirating capillary 6.1803.040 Connected to the injection valve. The other end can be connected either to the sample degasser or guided directly out of the instrument

9 Sample aspirating capillary 6.1803.040 Connected to the injection valve. Can optionally be connected to the sample degasser.

11 PEEK pressure screws, short 6.2744.070

For connecting the capillaries to the guard or separation column, respectively.

13 Conductivity detectors

8 PTFE connection capillary

Section of the 6.1803.040 PTFE capillary. Connects the sample degasser with the Sample Processor (required only if the sample degasser is used)

10 PTFE connection capillary

Section of the 6.1803.040 PTFE capillary. Connects the sample degasser with the Sample Processor (required only if the sample degasser is used)

12 Separation columns

You will find detailed descriptions of the individual installation steps in the following chapters.

3.4 Setting up the instrument

3.4.1 Packaging

The instrument is supplied in protective packaging together with the separately packed accessories. Keep this packaging, as only this ensures safe transportation of the instrument.

3.4.2 Checks

Immediately after receipt, check whether the shipment has arrived complete and without damage by comparing it with the delivery note.

3.4.3 Location

The instrument has been developed for operation indoors and may not be used in explosive environments.

Place the instrument in a location of the laboratory which is suitable for operation and free of vibrations and which provides protection against corrosive atmosphere and contamination by chemicals.

The instrument should be protected against excessive temperature fluctuations and direct sunlight.

3.5 Capillary connections in the IC system

This chapter contains general information concerning the capillary connections in the IC instruments and systems.

Generally speaking, capillary connections between two components of an IC system are made up of one connection capillary and two pressure screws with which the capillary is connected to the respective components.

Pressure screws

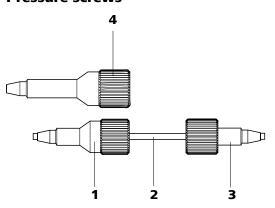


Figure 4 Connection of capillaries with pressure screws

- **1 PEEK pressure screw (6.2744.014)** Use on the injection valve.
- **3** PEEK pressure screw, short (6.2744.070)

For use on the high pressure pump, the purge valve, the inline filter, the pulsation absorber, the guard column and the separation column.

2 Connection capillary

4 PEEK pressure screw, long (6.2744.090)

Use on special components. Is not used on all instruments.



NOTICE

In order to keep the dead volume as low as possible, capillary connections should generally be as short as possible.



NOTICE

For an improved overview, capillary and tubing connections can be bundled with the 6.1815.010 spiral band.

Connection capillaries

PEEK capillaries and PTFE capillaries are used in the IC system.

PEEK capillaries (polyetheretherketone) PEEK capillaries are temperature-resistant up to 100°C, stable under pressure up to 400 bar, flexible, chemically inert and exhibit an extremely smooth surface. They can be readily cut down to the desired length with the 6.2621.080 capillary cutter.

Usage:

- PEEK capillaries (6.1831.010) with an internal diameter of 0.25 mm for the entire high-pressure section.
- PEEK capillaries (6.1831.030) with an internal diameter of 0.75 mm for sample handling in the ultratrace range.



CAUTION

For the capillary connections between the injection valve and detector, PEEK capillaries with an internal diameter of 0.25 mm must be used. These are already connected to a newly delivered instrument.

PTFE capillaries (polytetrafluoroethylene)

PTFE capillaries are transparent and enable visual tracing of the liquids to be pumped. They are chemically inert, flexible and temperature-resistant up to 80°C.

Usage:

PTFE capillaries (6.1803.0x0) are used for the low-pressure section.

- PTFE capillaries with internal diameter of 0.5 mm for sample handling.
- PTFE capillaries with internal diameter of 0.97 mm for sample handling as well as for rinsing solutions (they do not have to be in the scope of delivery of the instrument).

Capillary connections

In order to achieve optimum analysis results, capillary connections in an IC system must be absolutely tight and free of dead volume. Dead volume occurs if two capillary ends connected to each other do not fit exactly, thus allowing liquid to escape. There are two possible reasons for this:

- The capillaries do not have exactly cut edges.
- The two capillary ends do not completely meet.

One prerequisite for dead volume free capillary connection is, that both capillary ends are cut exactly plane. Therefore we recommend only to cut PEEK capillaries with the capillary cutter (6.2621.080).

Creating dead volume free capillary connections

To create dead volume free capillary connections, proceed as follows:

- 1 Slide the pressure screw over the capillary. Ensure that the capillary protrudes 1–2 mm from the tip of the pressure screw.
- Plug the capillary all the way into the connection or coupling until the stop.
- Only then start turning the pressure screw, while keeping the capillary pressed in space.

Colored sleeves for PEEK capillaries

The enclosed set of varicolored sleeves for PEEK capillaries (6.2251.000) serves to easily differentiate the various flows of liquid in the system through color coding. Each capillary leading a given liquid (e. g. eluent) can be highlighted with sleeves of the same color.

To highlight a capillary, proceed as follows:

Slide a sleeve of a selected color over a capillary an move it to an easily visible position.

If the capillary heats up, the sleeve shrinks and adapts to the form of the capillary.

3.6 Rear of the instrument

3.6.1 Rollers and handle

In order to make transport easier, the instrument is equipped with rollers and a handle.

3.6 Rear of the instrument

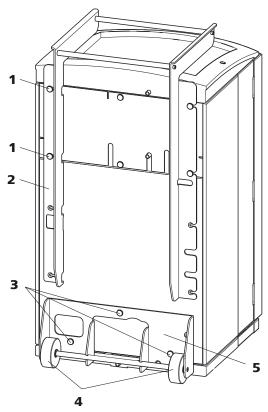


Figure 5 Rollers and handle

- **1 Knurled screws**For fastening the handle (5-2) and the rear panel of the detector chamber.
- **3** Knurled screws
 For fastening the roller holder (5-**5**).
- 5 Roller holder

- 2 Handle
- 4 Rollers

Removing handle

1 Loosen knurled screws (5-**1**) and remove handle (5-**2**).

Removing rollers

Proceed as follows to remove the rollers:

- **1** Remove knurled screws (5-**3**).
- **2** Remove roller holder (5-**5**).

Mounting handle as MPak holder



NOTICE

When extended, the handle (6-2) can also be used for hanging up MPaks (eluent bags).

1 Move handle (6-2) upwards and screw in the knurled screws (6-1) again.

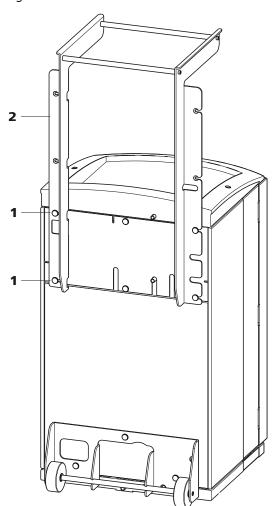


Figure 6 Handle as MPak holder

1 Knurled screws

For fastening the handle (6-2) and the rear panel of the detector chamber.

2 Handle

Extended. As holder for MPaks (eluent bag).

3.6 Rear of the instrument

3.6.2 Positioning and connecting the detector

The instrument is delivered without detector. For information on how to place and connect the detector, please refer to the manual of the detector.

3.6.3 Transport locking screws

To avoid damage to the high pressure pump and the vacuum pump during transport, the pumps are secured with transport locking screws.

Remove these transport locking screws before the initial start-up.

Removing transport locking screws

1 Remove all of the transport locking screws with the 4 mm hexagon key (6.2621.030) and keep them in a safe place.



WARNING

In order to avoid damage to the pumps, the transport locking screws must be remounted each time the instrument undergoes major transport.

3.6.4 Leak sensor

The leak sensor detects escaping liquid which collects in the base tray of the instrument.

To activate the leak sensor, the leak sensor connector plug (7-2) must be connected, the instrument switched on and the leak sensor switched to active in the software.

Connecting the leak sensor

Plug the leak sensor connector plug (7-2) into the leak sensor connection socket (7-1) on the rear of the instrument (see figure 7, page 21).

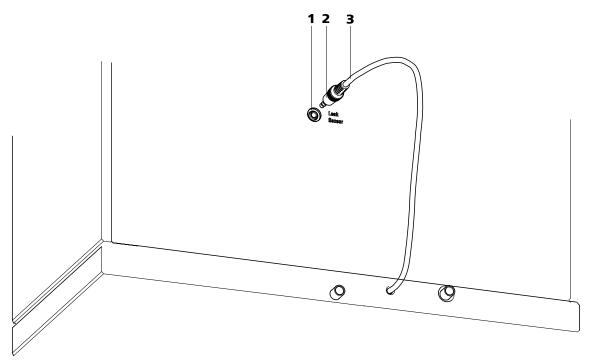


Figure 7 Connection of the leak sensor on the rear of the instrument

- 1 Leak sensor connection socket Is labeled with "Leak Sensor".
- 3 Leak sensor connection cable
 Is firmly mounted on the rear of the instrument.

2 Leak sensor connector plug

3.6.5 Drainage tubings

Liquid that escapes in the flask holder or in the detector chamber is directed through the drainage tubings into the base tray, past the leak sensor into the waste container. This ensures that any leaks in the system will be detected by the leak sensor.

3.6 Rear of the instrument

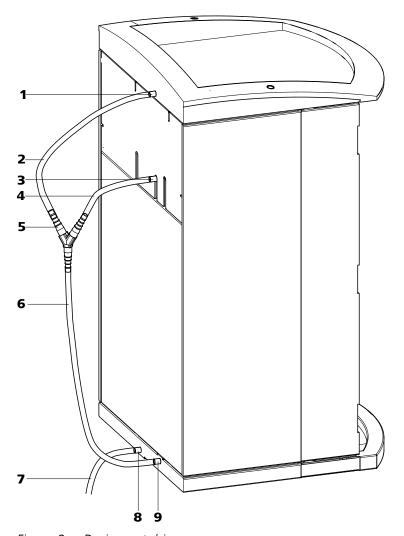


Figure 8 Drainage tubings

1 Drainage tubing connector

For draining escaped liquid from the flask holder.

3 Drainage tubing connector

For draining escaped liquid from the detector chamber.

5 Y connector 6.1807.010

For connecting the two drainage tubings (8-2) and (8-4).

2 Drainage tubing

Section of the 6.1816.020 silicon tubing. For draining escaped liquid from the flask holder.

4 Drainage tubing

Section of the 6.1816.020 silicon tubing. For draining escaped liquid from the detector chamber.

6 Drainage tubing

Section of the 6.1816.020 silicon tubing. Guides escaped liquid to the leak sensor.

' Drainage tubing

Section of the 6.1816.020 silicon tubing. Guides escaped liquid into a waste container

9 Drainage tubing connector

For supplying escaped liquid through the connected drainage tubing to the leak sensor.

8 Drainage tubing connector

For draining the escaped liquid from the base tray through the connected drainage tubing.

Proceed as follows to install the drainage tubings:

Installing drainage tubings

- Plug drainage tubing (8-2) into the drainage tubing connector (8-1) of the flask holder and shorten to required length.
- Plug drainage tubing (8-4) into the drainage tubing connector (8-3) of the detector chamber and shorten to required length.
- Connect drainage tubing (8-2) from the flask holder and drainage tubing (8-4) from the detector chamber to the Y connector (8-5).
- 4 Connect drainage tubing (8-6) to the Y connector (8-5), shorten to required length and plug the other end of the drainage tubing into the drainage tubing connector (8-9) of the base tray.
- Plug drainage tubing (8-7) into the drainage tubing connector (8-8) of the base tray and guide the other end into a waste container.

3.7 Capillary and cable feed-throughs

Several openings have been integrated for feeding through capillaries and cables. They are located on the doors (see figure 9, page 24), on the rear panel or below the flask holder or above the base tray (see Figure 10, Page 25).

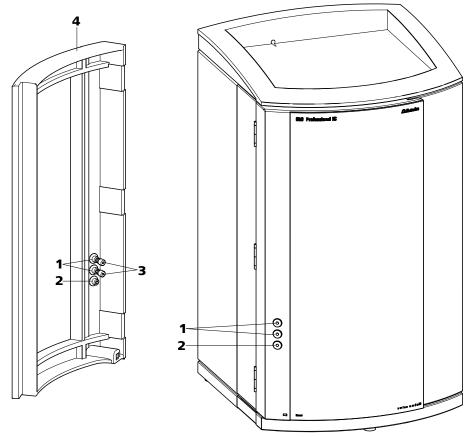


Figure 9 Capillary feed-throughs on the door

- **1 Luer connectors**For connecting a 6.2816.020 syringe. For manual sample feeding.
- **3** PEEK pressure screws, short 6.2744.070
- 2 Capillary feed-through
- 4 Door

Do not feed capillaries through the Luer connectors (9-1). The capillaries are fastened with PEEK pressure screws (9-3) from inside to the Luer connector. From outside, liquid can be aspirated or injected with a syringe.

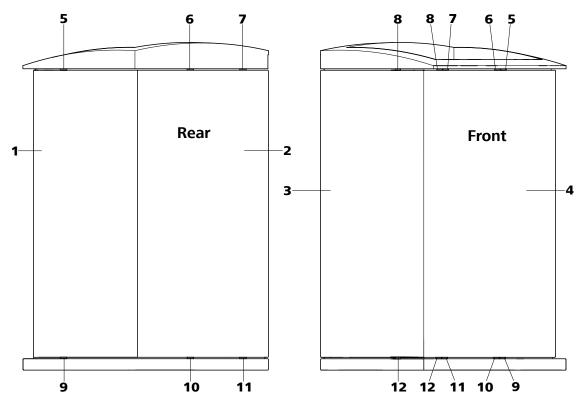


Figure 10 Capillary feed-throughs base tray/flask holder

1	Side panel (right) Right panel.	2	Rear of the instrument
3	Side panel (left) Left panel.	4	Front of the instrument
5	Capillary feed-through Upper. From front to right.	6	Capillary feed-through Upper. From front to back.
7	Capillary feed-through Upper. From front to back.	8	Capillary feed-through Upper. From front to left.
9	Capillary feed-through Lower. From front to right.	10	Capillary feed-through Lower. From front to back.
11	Capillary feed-through Lower. From front to back.	12	Capillary feed-through Lower. From front to left.

3.8 Eluent

3.8 Eluent

3.8.1 Connecting the eluent bottle

The eluent is aspirated out of the eluent bottle via the eluent aspiration tubing (11-1).

The eluent aspiration tubing is connected to the eluent degasser (see chapter 3.9, page 30). The tubing must be threaded through a suitable capillary feed-through (see chapter 3.7, page 23) of the instrument before the other end can be equipped.

The following accessories are needed for equipping the eluent aspiration tubing:

- 6.1602.160 eluent bottle cap GL 45
- 6.2744.210 tubing adapter for aspiration filter
- 6.2821.090 aspiration filter

Proceed as follows for equipping the eluent aspiration tubing:

Assembling eluent aspiration tubing

1 Guide the free end of the eluent aspiration tubing (11-1) out of the instrument through a suitable capillary feed-through.

2 Installing the eluent bottle cap (6.1602.160)

- Slide tubing nipple (11-2) and O-ring (11-3) onto the eluent aspiration tubing (11-1).
- Push eluent aspiration tubing (11-**1**) through the bottle cap (11-**4**) and screw tight.

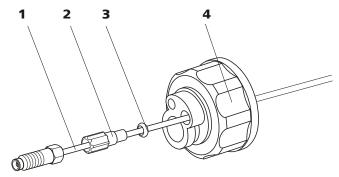


Figure 11 Installing the eluent bottle cap

- 1 Eluent aspiration tubing (6.1834.080)
- **Tubing nipple** From accessory set (6.1602.160).

3 O-ring

From accessory set (6.1602.160).

4 Bottle cap

From accessory set (6.1602.160).

3 Mounting the aspiration filter

• Insert the filter holder (12-1) into the aspiration filter (12-2) and screw it tightly.

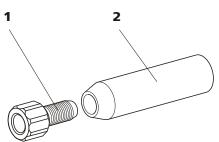


Figure 12 Mounting the aspiration filter

1 Filter holder

From accessory set (6.2744.210).

2 Aspiration filter (6.2821.090)

3.8 Eluent

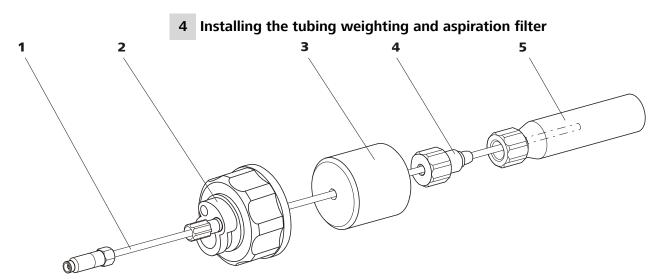


Figure 13 Installing the tubing weighting and aspiration filter

(6.2744.210).

1	Eluent aspiration tubing (6.1834.080)	2	2 Eluent bottle cap (6.1602.160)
3	Tubing weighting From accessory set (6.2744.210).	4	Clamping screw From accessory set (6.2744.210).
5	Aspiration filter (6.2821.090) With filter holder from accessory set		

- Slide the tubing weighting (13-3) onto the eluent aspiration tubing (13-1).
- Slide the clamping screw (13-4) onto the eluent aspiration tubing (13-1).
- Insert eluent aspiration tubing (13-1) into the aspiration filter (13-5). The end of the tubing should reach approximately to the center of the aspiration filter.
- Screw the clamping screw (13-4) to the filter holder (12-1).

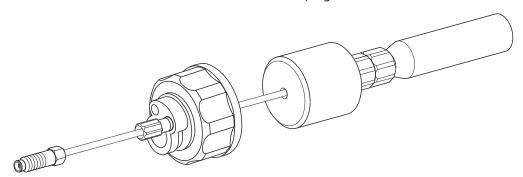


Figure 14 Eluent aspiration tubing fully equipped

5 Mounting the eluent aspiration tubing to the eluent bottle

• Insert the eluent aspiration tubing into the eluent bottle (15-10).

■ Tighten the fully equipped bottle cap to the eluent bottle (15-10). The aspiration filter (15-6) must rest on the base of the eluent bottle.

• Close the remaining small opening on the bottle cap with a threaded stopper (15-14) from the accessory set.

6 Mounting the adsorber tube



NOTICE

If alkaline eluents or eluents with lower buffer capacity are used, the eluent bottle must be equipped with an adsorber tube filled with CO_2 adsorber (15-**4**).

- First, place a piece of cotton (15-3), then the CO₂ adsorber (15-4) in the large opening of the adsorber tube (15-2) and close with the plastic cover.
- Fasten the adsorber tube (15-2) on the bottle cap (15-11) using the SGJ clip (15-12).

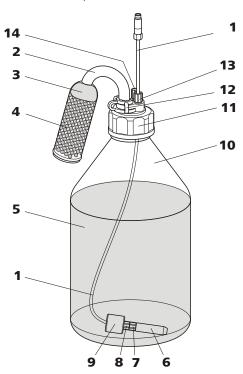


Figure 15 Eluent bottle – connected

1 Eluent aspiration tubing (6.1834.080) For aspirating the eluent. Pre-installed.

2 Adsorber tube (6.1609.000)

3.9 Eluent degasser

3	Cotton	4	CO ₂ adsorber Adsorbs CO ₂ from the air (e.g. Merck soda lime with indicator, no. 6839.1000).
5	Eluent	6	Aspiration filter (6.2821.090)
7	Filter holder From accessory set (6.2744.210).	8	Clamping screw From accessory set (6.2744.210).
9	Tubing weighting From accessory set (6.2744.210).	10	Eluent bottle (6.1608.070)
11	Bottle cap (6.1602.160)	12	SGJ clip (6.2023.020)
13	Tubing nipple	14	Threaded stopper

3.9 Eluent degasser

Gas bubbles in the eluent lead to an unstable baseline, as high pressure pumps can transport liquids, but not gases. This is why the eluent must be degassed before it reaches the high-pressure pump.

The eluent degasser removes gas bubbles and dissolved gases from the eluent. For degassing, the eluent flows into a vacuum chamber through a special fluoropolymer capillary.



NOTICE

The eluent degasser is already installed in the newly delivered instrument. The following installation instructions must only be followed, if the connections to the degasser had to be disconnected for maintenance.

Connecting the eluent degasser

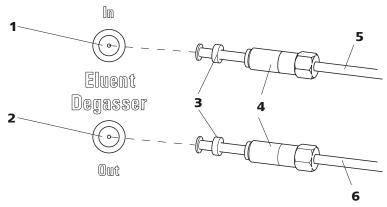


Figure 16 Eluent degasser

- 1 Eluent degasser input
- **Tubing flare**With tubing nipple.
- **5 Eluent aspiration tubing (6.1834.080)**For aspirating the eluent. The clamping screw (16-**4**) is firmly mounted.
- 2 Eluent degasser output
- 4 Clamping screw
- **Connection tubing (6.1834.090)**Connection from the eluent degasser to the high-pressure pump (see chapter 3.10, page 32). The clamping screw (16-**4**) is firmly mounted.

1



CAUTION

The clamping screws (16-4) must be tightened carefully. Use the wrench (6.2621.050) to do this.

- Insert the eluent aspiration tubing (16-**5**) into the eluent degasser input (16-**1**).
- Carefully tighten the clamping screw (16-**4**).
- Insert connection tubing (16-6) (the end with the longer clamping screw (16-4)) into the eluent degasser output (16-2).
 - Carefully tighten the clamping screw (16-**4**).
 - Connect the other end of the connection capillary (16-6) (with the shorter clamping screw) to the high-pressure pump (17-9) (see "Connecting inlet to the high pressure pump", page 33).

3.10 High pressure pump

3.10 High pressure pump

The intelligent and low pulsation high pressure pump pumps the eluent through the system. It is equipped with a chip on which its technical specifications and "life history" (operating hours, service data, ...) are saved.

The purge valve is used for deaerating (see chapter 3.10.2, page 34) the high pressure pump.

3.10.1 Capillary connections high pressure pump/purge valve



NOTICE

All of the capillary connections of the high pressure pump and the purge valve are already installed in the newly delivered instrument.

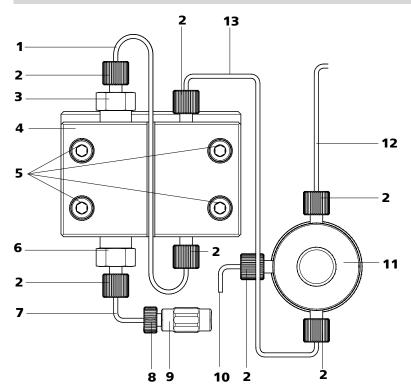


Figure 17 Capillary connections high pressure pump/purge valve

- **1** Connection capillary
 PEEK capillary, connects main piston and auxiliary piston.
- **3** Outlet valve holder

- 2 PEEK pressure screw, short (6.2744.070)
- 4 Pump head (6.2824.110)

5 Fastening screws

For fastening the pump head.

7 Pump head inlet capillary

PEEK capillary at the input of the pump head

9 Coupling

For the connection of the eluent path at the input of the high pressure pump. Can be ordered together with the pressure screw (17-8) under the number (6.2744.230).

11 Purge valve

For deaerating the high pressure pump. With rotary knob in the center and pressure sensor.

13 Connection capillary

Connects the output of the pump head with the purge valve.

6 Inlet valve holder

8 Pressure screw

For connecting a PEEK capillary to the coupling (17-**9**).

10 Deaerating capillary

For aspirating the eluent when deaerating the high pressure pump (see chapter 3.10.2, page 34).

12 Connection capillary

For connecting the inline filter (see chapter 3.11, page 36).



NOTICE

The eluent aspiration capillary is already installed in the newly delivered instrument. The following installation instructions need **not** be carried out at the time of initial installation.

Connecting inlet to the high pressure pump

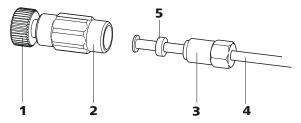


Figure 18 High pressure pump – Connect inlet

1 Pressure screw

For connecting the coupling (18-2) to the pump head inlet capillary (17-7).

Can be ordered together with the coupling under the number (6.2744.230).

2 Coupling (6.2744.230)

For connecting the eluent connection capillary (18-**4**) to the input of the high pressure pump.

3.10 High pressure pump

3 Clamping screw

4 Eluent aspiration tubing

Eluent aspiration tubing (6.1834.080) or (6.1834.090).

5 Backup ring

1 Connecting coupling

Fasten the coupling (18-2) with a pressure screw (18-1) on the pump head inlet capillary (17-7).

2 Connecting eluent aspiration tubing



CAUTION

The clamping screws must be tightened carefully. To tighten, grip the coupling (18-2) with the key (6.2739.000) and grip the clamping screw (18-3) with the wrench (6.2621.050).

- Plug the eluent aspiration tubing (18-4) into the coupling (18-2).
- Tighten the clamping screw (18-3).

3.10.2 Deaerating the high pressure pump

The high pressure pump will only operate perfectly if the pump head contains no more air bubbles. Therefore it must be deaerated during initial start-up and after every change of eluent.



CAUTION

The high pressure pump must **not** be deaerated before the initial startup .

Deaerate the high pressure pump as follows (see figure 19, page 35):

Deaerate the high pressure pump

The instrument must be connected to the PC and switched on to deaerate the high pressure pump.

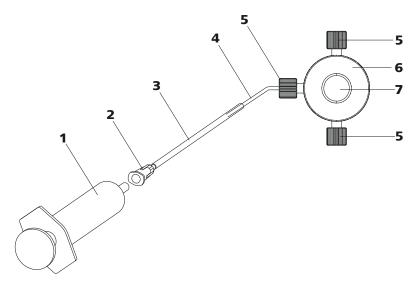


Figure 19 Deaerate the high pressure pump

1	Syringe 10 mL (6.2816.020) For aspirating the eluent.	2	Luer connector Part of the purging needle (6.2816.040)
3	Purging needle (6.2816.040)	4	Deaerating capillary
5	PEEK pressure screws, short (6.2744.070)	6	Purge valve
7	Purge valve rotary knob		

1 Connecting the purging needle

• Push the end of the purging needle (19-3) over the end of the deaerating capillary (19-4) on the purge valve.

2 Connecting the syringe

• Insert syringe (19-**1**) in the Luer connector (19-**2**) of the purging needle (see figure 19, page 35).

3 Opening purge valve

• Open the rotary knob (19-**7**) by approx. ½ rotation counterclockwise.

4 Setting the flow rate

- Start MagIC Net[™] (if not yet started).
- Ensure that the eluent aspiration tubing is immersed sufficiently in the eluent.
- Let the high pressure pump run.

3.11 Inline filter

5 Aspirating eluent

• Aspirate with the syringe (19-1) until bubble-free eluent flows into the syringe.

6 Completing deaerating

- Switch off high pressure pump.
- Close rotary knob (19-**7**).
- Remove syringe (19-1) from the Luer connector (19-2).
- Pull the purging needle (19-3) out of the deaerating capillary (19-4).

3.11 Inline filter

Between the purge valve and the pulsation absorber the inline filter (6.2821.120) is installed as protection against particles.

Inline filters protect the separation column securely against possible contamination from the eluent. Inline filters can however also just as well be used for the purpose of protecting the suppressor against contaminations in the regeneration or rinsing solutions. The filter platelets with a pore size of 2 μm can be replaced quickly and easily. They remove particles like e. g. bacteria and algae from the solutions.



NOTICE

The inline filter is already installed in the newly delivered instrument. The following installation instructions need **not** be carried out at the time of initial installation.

Installing the inline filter



CAUTION

Observe the flow direction marked on the filter housing for the connection of the inline filter.

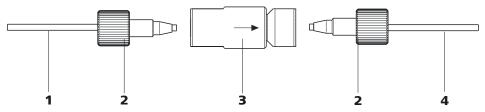


Figure 20 Connecting the inline filter

1 Connection capillary

Connects the purge valve with the inline filter

3 Inline filter (6.2821.120)
Protects against particles.

- 2 PEEK pressure screws, short (6.2744.070)
- 4 Connection capillary

Connects the inline filter with the pulsation absorber.

- 1 Screw on the connection capillary running from the purge valve to the input side of the inline filter using a pressure screw (6.2744.070).
- 2 Screw on the connection capillary running to the pulsation absorber to the output side of the inline filter using a pressure screw (6.2744.070).

3.12 Pulsation absorber



NOTICE

The pulsation absorber is already installed in the newly delivered instrument.



CAUTION

The pulsation absorber is maintenance-free and may not be opened.

The pulsation absorber protects the separation column from damage caused by pressure fluctuations when switching the injection valve, and reduces interfering pulsations during highly sensitive measurements. In order to ensure these functionalities, it must be connected between the high pressure pump (see chapter 3.10, page 32) and injection valve (see chapter 3.14, page 40).

The pulsation absorber can be operated in both directions.

3.13 Sample degasser

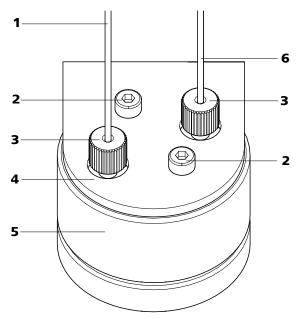


Figure 21 Pulsation absorber – Connection

1 Connection capillary
Connection to the inline filter.

2 Fastening screws

3 PEEK pressure screws, short
(6.2744.070)

4 Holder for pulsation absorber

6 Connection capillary
Connection to the injection valve.

3.13 Sample degasser

The sample degasser removes gas bubbles and dissolved gases from the sample. For degassing, the sample flows into a vacuum chamber through a special fluoropolymer capillary.

Gas bubbles in the sample lead to poor reproducibility, as the amount of sample in the sample loop would not always be the same. This is why samples (containing gas) should be degassed prior to injection. The sample is drawn in through a degassing chamber prior to injection and any gas bubbles are automatically removed.



NOTICE

The rinsing time increases by at least 2 minutes when the sample degasser is used.

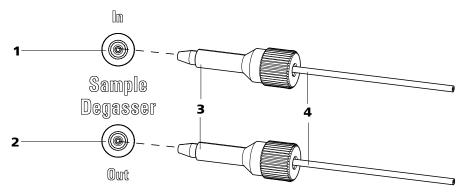


Figure 22 Sample degasser

1 Sample de	egasser input	2	Sample degasser output
3 PEEK pres (6.2744.09	sure screw, long 90)	4	Connection capillaries (6.1803.040)

Connecting the sample degasser

- Remove and keep the threaded stoppers (6.2744.220) from the inlet and outlet of the sample degasser.
- 2 Connect the end of the sample aspiration capillary (6.1803.040) connected to the injection valve to the output of the sample degasser (22-2), using a long PEEK pressure screw (22-3).
- **3** Connect the connection capillary (6.1803.040) to the input of the sample degasser (22-1), using a long PEEK pressure screw (22-3).
- **4** Guide the other end of the connection capillary out of the instrument through a capillary feed-through and connect it with the Sample Processor, if applicable.



CAUTION

If the sample degasser is not used, the input and output **must** be sealed with the threaded stoppers (6.2744.220).

3.14 Injection valve

3.14 Injection valve

The injection valve connects the eluent and sample path. Through rapid and precise valve switchover a precise amount of sample solution defined by the size of the sample loop is injected and rinsed with eluent onto the separation column.

3.14.1 Connecting the injection valve

The injection valve has six connectors: two for the sample path (connectors 1 and 2), two for the eluent path (connectors 4 and 5) and two for the sample loop (connectors 3 and 6).



NOTICE

The capillaries of the eluent path and the sample path and the sample loop are already installed in the newly delivered instrument.

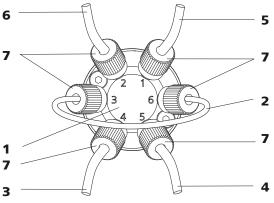


Figure 23 Injection valve – connected

1 Injection valve

3 Connection capillary

Connected to connector 4. Carries eluent to the injection valve.

5 Connection capillary

Connected to connector 1. Carries sample to the injection valve.

7 PEEK pressure screw 6.2744.010

2 Sample loop

Connected to connectors 3 and 6.

4 Connection capillary (column inlet capillary)

Connected to connector 5. Carries eluent to the separation column.

6 Connection capillary

Connected to connector 2. Carries sample to the waste container.

Replacing the sample loop

The sample loop can be replaced, depending on requirements. For additional information concerning selection of the appropriate sample loop, see Chapter 3.14.3, page 42.



NOTICE

Use only 6.2744.010 PEEK pressure screws for connecting capillaries and sample loop to the injection valve.

1 Removing existing sample loop

- Loosen 6.2744.010 pressure screws at connector 3 and connector
 6.
- Remove sample loop.

2 Mounting new sample loop

- Fasten one end of the sample loop (23-2) with a 6.2744.010 PEEK pressure screw (23-7) to connector 3.
- Fasten the other end of the sample loop (23-2) with a second 6.2744.010 PEEK pressure screw (23-7) to connector 6.

3.14.2 Mode of operation of the injection valve

The injection valve (see figure 24, page 41) can adopt two valve positions - **FILL** and **INJECT**. Switching back and forth between the two valve positions determines whether the sample path or the eluent path is guided through the sample loop. The following figure provides a schematic display of the flow paths of the two valve positions.

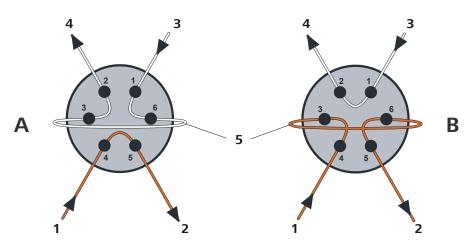


Figure 24 Injection valve – Positions

A Position FILL

B Position INJECT

3.15 Column thermostat

1 Eluent input

Capillary coming from the high pressure pump.

3 Sample input

Sample aspirating capillary.

5 Sample loop

2 Eluent output

Capillary to the column.

4 Sample output

Capillary to waste container.

Position A

In the position **FILL**, the sample solution flows through the sample loop to the waste container. The eluent flows directly to the separation column at the same time.

Position B

In the position **INJECT**, the eluent flows through the sample loop to the separation column. If sample solution is to be found in the sample loop at the time of the valve switchover, then this will be conveyed along with the eluent, thus making its way to the separation column. The flow in the sample path is either stopped or the sample flows directly to the waste container.

3.14.3 Selecting the sample loop

The amount of sample solution injected depends on the volume of the sample loop. The choice is made on the basis of the application. The following sample loops are normally used:

Cation determination 10 μ L

Anion determination with suppression 20 μ L

Anion determination without suppression 100 μ L

3.15 Column thermostat

The column thermostat controls the temperature of the column and eluent channel and thus ensures stable measuring conditions. It provides space for 2 separation columns.

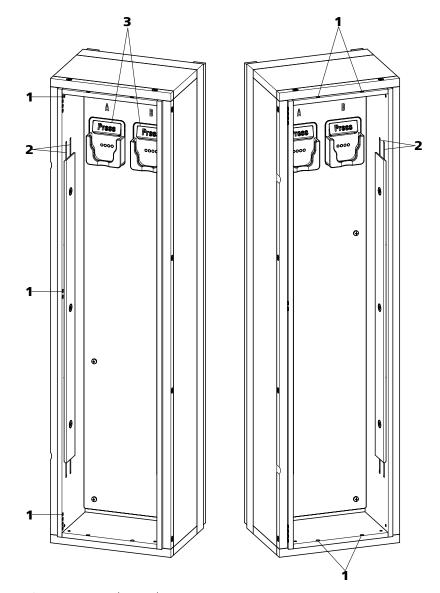


Figure 25 Column thermostat

1 Capillary feed-throughs

For guiding the capillaries in and out.

2 Capillary recesses

For controlling the temperature of the eluent.

Preheating capillary already pre-installed.

3 Column holder

For fastening the columns.

With column recognition.

The column thermostat contains two column holders (25-**3**) equipped with chip recognition. The separation columns can be clicked into the column holder with the chip.

3.15 Column thermostat



NOTICE

The column input capillary is already threaded into the capillary recesses of the column thermostat at the time the instrument is first delivered. The following installation instructions need **not** be carried out at the time of initial installation.

Threading capillaries

- 1 Introduce column input capillary into the column thermostat via a suitable capillary feed-through (25-1).
- 2 Slide the column input capillary from below into the outer of the two capillary recesses (25-2). Slide it through under the holder plate until it emerges again at the top.
- **3** Carefully bend the column input capillary downward and slide from from above to below through the inner capillary recess until it emerges at the lower edge of the holder plate.

4



NOTICE

The columns (guard column and separation column) may only be installed after the initial start-up .

Before initial start-up:

Fasten the 6.2744.040 coupling with a 6.2744.010 pressure screw to the end of the column input capillary.

• After initial start-up:

Fasten the guard column (if used) or the separation column with a 6.2744.010 pressure screw to the end of the column input capillary.

3.16 Conductivity detector

The conductivity detector continuously measures the conductivity of the liquid passing through and indicates these signals in digital form (DSP – Digital Signal Processing). The conductivity detector exhibits outstanding thermal stability and thus guarantees reproducible measuring conditions.

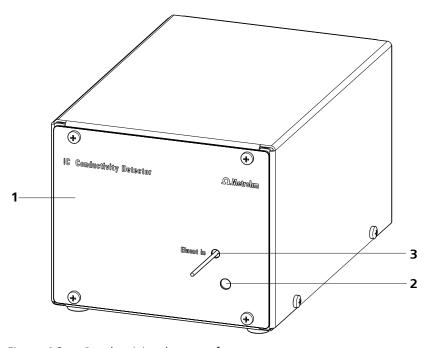


Figure 26 Conductivity detector front

1 IC detector 1.850.9010

2 Opening for temperature sensor

3 Detector input capillary Permanently installed.

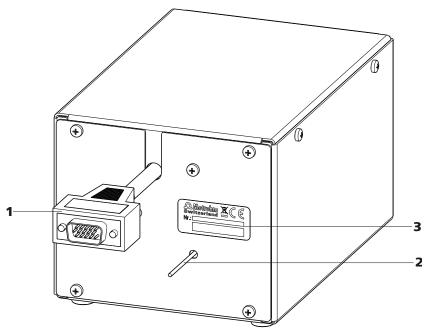


Figure 27 Conductivity detector rear

- **1 Detector cable**With mounted plug.
- **Type plate**With serial number.

2 Detector output capillary Permanently installed.



NOTICE

In order to prevent unnecessary peak widening after separation, the connection between the outlet of the separation column and the inlet in the detector should be kept as short as possible.

Connect the detector input capillary to the separation column

- 1 Connecting detector input
 - Fasten the detector input capillary (28-1) on the output of the column (28-3) using a 6.2744.070 pressure screw (28-2).

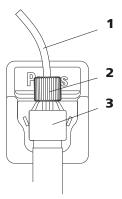


Figure 28 Connection detector – separation column

1 Detector input capillary

2 PEEK pressure screw, short 6.2744.070

3 Separation column

3.17 Connecting the instrument to a computer



NOTICE

If the instrument is connected to the computer, then it must be switched off.

Accessories

For this step, you need the following accessories:

■ USB connecting cable (6.2151.020)

Connecting the USB cable

- 1 Insert the USB cable into the computer connection socket on the rear of the instrument labeled *PC*.
- 2 Insert the other end into a USB port on the computer.

3.18 Connecting the instrument to the power grid



WARNING

Electric shock from electrical potential

Risk of injury by touching live components or through moisture on live parts.

- Never open the housing of the instrument while the power cord is still connected.
- Protect live parts (e.g. power supply unit, power cord, connection sockets) against moisture.
- Unplug the power plug immediately if you suspect that moisture has gotten inside the instrument.
- Only personnel who have been issued Metrohm qualifications may perform service and repair work on electrical and electronic parts.

Connecting the power cord

Accessories

Power cord with the following specifications:

- Length: max. 2 m
- Number of cores: 3, with protective conductor
- Instrument plug: IEC 60320 type C13
- Conductor cross-section 3x min. 0.75 mm² / 18 AWG
- Power plug:
 - according to customer requirement (6.2122.XX0)
 - min. 10 A



NOTICE

Do not use a not permitted power cord!

1 Plugging in the power cord

- Plug the power cord into the instrument's power socket.
- Connect the power cord to the power grid.

3.19 Guard column

The use of guard columns serves for protecting the separation columns and increasing their service life considerably. The guard columns available from Metrohm represent either actual guard columns or are so-called guard column cartridges which are used together with a cartridge holder. The installation of a guard column cartridge in the associated holder is described in the leaflet of the guard columns.



NOTICE

Information regarding which guard column is suitable for your separation column can be found in the **Metrohm IC Column Program** (which is available from your Metrohm agent), the leaflet provided along with your separation column, the product information on the separation column at http://www.metrohm.com (product area Ion Chromatography), or obtained directly from your agent.



CAUTION

New guard columns are filled with solution and are sealed on both sides with stoppers or caps, respectively. Before using the guard column, you need to ensure that this solution is miscible with the eluents used (observe manufacturer's data).



NOTICE

The guard column may only be installed after the **initial start-up** (see chapter 4.1, page 53) of the instrument. Until then, use the coupling (6.2744.040) instead of the guard and separation column.



NOTICE

Metrohm recommends, always to work with guard columns. These protect the separation column and can be exchanged regularly.

3.20 Separation column

Connecting and rinsing the guard column

1 Connecting the guard column



CAUTION

When inserting the guard column, always ensure that it is inserted correctly corresponding to the flow direction (if indicated).

- Remove sealing caps and/or stoppers from the guard column.
- Fasten the input of the guard column to the column inlet capillary using a short PEEK pressure screw (6.2744.070).
- In case the guard column is mounted on the separation column with a connecting capillary, connect the included connection capillary, which is included with the guard column, to the output of the guard column using the PEEK pressure screw, which is .also included.

2 Rinsing the guard column

- Place beaker under the outlet capillary of the guard column.
- Set the flow rate of the high pressure pump according to the data given in the leaflet of the separation column.
- Start the high pressure pump and rinse the guard column approx.
 5 minutes with eluent.
- Switch off the high pressure pump again.

3.20 Separation column

The intelligent separation column (iColumn) is the heart of the ion chromatographic analysis. It separates the different components corresponding to their interactions with the column. Metrohm separation columns are equipped with a chip on which their technical specifications and their history (first use / setting up, operating hours, injections, ...) are saved.



NOTICE

Information regarding which separation column is suitable for your application can be found in the **Metrohm IC Column Program**, the product information for your separation column at http://www.metrohm.com in the product area ion chromatography, or obtained directly from your agent.



CAUTION

New separation columns are filled with solution and are sealed on both sides with stoppers. Before using the column, you need to ensure that this solution is miscible with the eluents used (observe manufacturer's data).

You can find the separation columns and guard columns currently available from Metrohm in the Metrohm IC Column Program, or in the Internet at http://www.metrohm.com in the product area Ion Chromatography. A test chromatogram and an leaflet are provided along with each column. You can request detailed information on special IC applications in the corresponding "Application Bulletins" or "Application Notes", available in the Internet at http://www.metrohm.com in the Applications area or via the Metrohm agent responsible free of charge.



NOTICE

The separation column may only be installed after the **initial start-up** (see chapter 4.1, page 53) of the instrument. Until then, use the coupling (6.2744.040) instead of the guard and separation column.

Connecting and rinsing the separation column

1 Connect the separation column



CAUTION

When inserting the columns, always ensure that these are correctly inserted corresponding to the flow direction indicated.

- Remove stoppers from the separation column.
- Attach the guard column to the input of the separation column.
 OR

Connect the input of the separation column to the outlet capillary of the guard column, using the PEEK pressure screw (6.2744.070) included.

OR

If no guard column is used (not recommended), connect the column input capillary to the input of the separation column, using a PEEK pressure screw (6.2744.070).

3.20 Separation column

2 Rinsing the separation column

- Place beaker under the outlet end of the separation column.
- Set the flow rate of the high pressure pump according to the data given in the leaflet of the separation column.
- Start the high pressure pump and rinse the separation column approx. 10 minutes with eluent.
- Switch off the high pressure pump again.

3 Mounting the separation column

- Fasten the column output capillary to the output of the separation column using a PEEK pressure screw (6.2744.070).
- Hang separation column with chip into the column holder.



NOTICE

The iColumns are equipped with a chip on which their operating data is saved. The chip has to be hooked into the chip holder provided for this so that the column recognition can function.

4 Start-up

4 Start-up

The chapter *Start-up* is divided into 2 sections:

Initial start-up The initial start-up is carried out during the ini-

tial installation.

Conditioning Conditioning is carried out as a final installation

step and each time after the system is started.

4.1 Initial start-up

The initial start-up is carried out during the initial installation. The entire system is rinsed before guard column and separation column are installed.



CAUTION

The separation column and guard column may not be installed for the initial start-up.

Make sure that the coupling (6.2744.040) is being used instead of the columns.

Perform the following steps during the initial start-up:

1 Preparing the software

- Start the PC program **MagIC Net**TM.
- Open the **Equilibration** tab in **MagIC Net**TM.
- Select (or create) a suitable method.

2 Preparing the instrument

- Ensure that the eluent aspiration tubing is immersed in the eluent and that there is enough eluent in the eluent bottle.
- Switch on the instrument.

3 Deaerate the high pressure pump

• Deaerate the high pressure pump(s) via the purge valve (see chapter 3.10.2, page 34).

4 Rinsing the instrument without columns

• Rinse the instrument (without columns) with eluent for 5 minutes.

4.2 Conditioning

The instrument is now ready for the installation of the columns.

4.2 Conditioning

After the installation and after switching on the instrument, the system must be conditioned with eluent until a stable baseline is reached.



NOTICE

After a change of eluent (see chapter 5.4.2.3, page 59), the conditioning time can lengthen considerably.

Conditioning the system

1 Preparing the software



CAUTION

Ensure that the flow set is not higher than the flow permissible for the corresponding column (see column leaflet and chip data set).

- Start the PC program MagIC Net™.
- Open the **Equilibration** tab in **MagIC Net**TM.
- Select (or create) a suitable method.

2 Preparing the instrument

- Ensure that the column is correctly mounted according to the flow direction indicated on the label (arrow must point in the direction of flow).
- Ensure that the eluent aspiration tubing is immersed in the eluent and that there is enough eluent in the eluent bottle.

3 Checking leak-tightness

- In **MagIC Net**[™], start the equilibration.
- Check all capillaries and their connections from the high pressure pump to the detector block for signs of liquid escaping. If eluent escapes anywhere, tighten the corresponding pressure screw or loosen the connection, check the end of the capillary, shorten it with a capillary cutter if necessary, and restore the connection.

4 Start-up

4 Conditioning the system

Rinse the system with eluent until the required stability of the base-line is attained (normally 30 minutes).

The instrument is now ready for measuring samples.

5.1 General notes

5 Operation and maintenance

5.1 General notes

5.1.1 Care



WARNING

The instrument housing must not be opened by untrained personnel.

The instrument requires appropriate care. Excess contamination of the instrument may result in functional disruptions and a reduction in the service life of the sturdy mechanics and electronics.



CAUTION

Although this is prevented to a great extent by design measures, the mains plug should be unplugged immediately if aggressive media has penetrated the inside of the instrument, so as to avoid serious damage to the instrument electronics. In such cases, the Metrohm Service must be informed.

On the rear of the instrument, the drainage tubings must be mounted and the leak sensor must be plugged in and activated as protection against escaping liquids.

Spillages of chemicals and solvents should be cleaned up immediately. In particular, the plug connections on the rear panel of the instrument (especially the mains plug) should be protected from contamination.

5.1.2 Maintenance by Metrohm Service

Maintenance of the instrument is best carried out as part of an annual service, which is performed by specialist personnel from Metrohm. If working frequently with caustic and corrosive chemicals, a shorter maintenance interval is recommended. The Metrohm service department offers every form of technical advice for maintenance and service of all Metrohm instruments.

5 Operation and maintenance

5.1.3 Operation



CAUTION

In order to avoid disturbing temperature influences, the entire system including the eluent bottle must be protected against direct sunlight.

5.1.4 Shutting down

If the instrument is not used for a longer period, the whole IC system (except the columns) must be rinsed salt free with methanol/ultrapure water (1:4), in order to prevent eluent salts from forming crystals which may cause subsequent damage.

Rinsing salt free the IC system

To rinse the system, proceed as follows:

- Remove the separation column from the eluent path. Connect the connection capillaries directly with each other using a coupling (6.2744.040).
- **2** Rinse the IC system with methanol/ultrapure water (1:4) for 15 minutes.

Rinse with eluent for at least 15 minutes at starting up again and before connecting the guard column and separation column.

5.2 Capillary connections

5.2.1 Operation

All connections between injection valve, separation column and detector must be as short as possible, have a low dead volume and be completely leak-tight. The PEEK capillary after the detector must be free of blockages. Only use PEEK capillaries with an internal diameter of 0.25 mm in the high-pressure section between the high pressure pump and the detector.

5.3 Door

5.3 Door



CAUTION

The door is made of PMMA (polymethylmetacrylate). It must never be cleaned with abrasive media or solvents.



CAUTION

Never use the door as a handle.

5.4 Eluent

5.4.1 Production

Chemicals used for manufacturing eluents must have a purity grade of at least "p.a.". Only ultrapure water (resistance > 18.2 M Ω *cm) may be used for dilution (this generally applies for reagents which are used in ion chromatography).

Newly manufactured eluents always have to be microfiltered (0.45 μm filter).

The composition of the eluent plays a critical role in chromatographic analysis:

Concentration An increase in the concentration generally leads

to shorter retention times and faster separation, but also to a higher background conductivity sig-

nal.

pH pH changes lead to shifts in dissociation equili-

bria and thus to changes in retention times.

Organic solvents Adding an organic solvent (e.g. methanol, ace-

tone or acetonitrile) to watery eluents generally

speeds up lipophilic ions.

5.4.2 Operation

5.4.2.1 Supply bottle

The supply bottle with the eluent must be connected as indicated in *chapter 3.8.1*, page 26. This is particularly important for eluents with volatile solvents (e.g. acetone).

Condensation in the eluent bottle must be avoided. Drop formation can change the concentration ratio in the eluent.

In the case of very sensitive measurements, we recommend that the eluent be stirred constantly with a magnetic stirrer (e. g. the 2.801.0010 with 6.2070.000).

5.4.2.2 Aspiration filter

To protect the IC system from foreign particles, we recommend aspirating the eluent via an aspiration filter (6.2821.090) (12-2). This aspiration filter must be replaced if it takes on a yellowish discoloration, and at least every 3 months.

5.4.2.3 Changing the eluent

Prevent any precipitations when changing the eluent. Immediately successive solutions must always be miscible. If the system has to be rinsed with organic solvents, use several solvents with rising or falling lipophilicity.

5.5 High pressure pump

5.5.1 Protection



CAUTION

The pump head is filled ex works with methanol/ultrapure water. It must be ensured that the eluent used is freely miscible with the solvent remaining in the pump head.

To protect the high pressure pump against **foreign particles**, we recommend that the eluent undergoes a **microfiltration** (filter 0.45 μ m) before being aspirated via a 6.2821.090 aspiration filter (see "Assembling eluent aspiration tubing", page 26).

Salt crystals between the piston and seal cause abrasion particles which can find their way into the eluent. These lead to contaminated valves, a rise in pressure and in extreme cases scratched pistons. It is therefore essential to ensure that **no precipitates** can occur (see chapter 5.4.2.3, page 59).

5.5 High pressure pump



CAUTION

In order to spare the pump seals, the pump should not be operated dry. Therefore ensure that the eluent supply is correctly connected and that there is enough eluent in the eluent bottle each time before switching on the pump.

5.5.2 Maintenance



CAUTION

Maintenance work on the high pressure pump may not be carried out unless the **instrument is switched off**.

Pump head maintenance

An unstable baseline (pulsation, flow fluctuations) is in many cases the result of contaminated valves (35-2), (35-3) or defective, leaking piston seals on the high pressure pump. Proceed as follows for cleaning contaminated valves and/or replacing worn parts such as pistons, piston seal and valves:

This maintenance work should be carried out at least once a year.

Removing the pump head

- 1 Switch off high pressure pump and wait until pressure is released.
- **2** Loosen the pressure screw on the inlet valve holder (17-**6**) and unscrew the pump head input capillary (17-**7**), the coupling (17-**9**), and the eluent aspiration tubing from the pump head.
 - In the process, eluent will spill. Hold the eluent aspiration tubing up high and allow the eluent to flow back into the eluent bottle.
- **3** Unscrew the pump head outlet capillary (17-**13**) from the pump head.
- 4 Remove pump head from the pump housing by loosening the 4 fastening screws (17-**5**) using the hexagon key (6.2621.030). The main piston is on the left (viewed from the front), and the auxiliary piston is on the right.

5 Operation and maintenance

Cleaning/replacing the zirconium oxide piston

Clean one piston after the other as follows:

1 Removing the piston cartridge from the pump head

Loosen the piston cartridge with a wrench and unscrew from the pump head by hand.

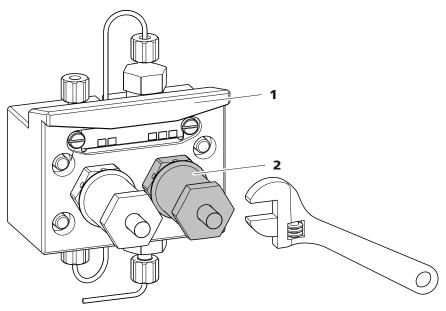


Figure 29 Pump head – removing the piston

1 Pump head

2 Piston cartridge

2 Dismantling the piston



CAUTION

On the inside of the piston cartridge there is a taut spring than can jump out of the piston cartridge if suddenly loosing tension.

When opening the piston cartridge, hold pressure towards the spring and unscrew carefully.

- Loosen the screw of the piston cartridge with a wrench and unscrew carefully by hand and by holding pressure towards the taut spring.
- Remove the zirconium oxide piston and lay on a tissue.
- Remove the spring retainer, spring and the inner plastic sleeve from the piston cartridge and lay by.

5.5 High pressure pump

• Remove the backup ring from the pump head and lay to the other parts.

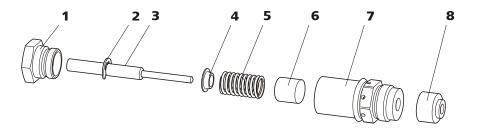


Figure 30 Components of the piston cartridge

1	Piston cartridge screw	2	Retaining washer
3	Zirconium oxide piston with piston shaft Order number: 6.2824.070	4	Spring retainer
5	Spring Order number: 6.2824.060	6	Inner plastic sleeve Protects from metallic abrasion.
7	Piston cartridge	8	Backup ring

3 Cleaning the components of the piston

- Clean zirconium oxide pistons contaminated by abrasion or deposits with fine abrasive cleaning powder, rinse particle free with ultrapure water and dry.
 Replace highly contaminated or scratched zirconium oxide pistons (spare part: 6.2824.070 zirconium oxide piston).
- Rinse the other parts of the piston and dry with a lint-free cloth.

4 Assembling the piston

- Insert the inner plastic sleeve, spring and spring retainer into the piston cartridge.
- Slide the zirconium oxide piston carefully into the piston cartridge until its tip emerges from the small opening of the piston cartridge.
- Attach screw and tighten by hand.

Replacing the piston seal

The special tool (6.2617.010) (see figure 31, page 63) is necessary to remove the piston seal from the pump head. It consists of two parts: a tip

5 Operation and maintenance

for removing the old piston seal and a sleeve for inserting the new piston seal.

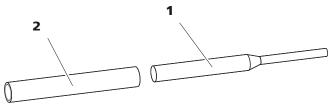


Figure 31 Tool for piston seal

1 Pin

Tip for removing the old piston seal.

2 Sleeve

Sleeve for inserting the new piston seal.



CAUTION

Screwing the special tool for the piston seal (6.2617.010) into the piston seal destroys this completely!

1 Removing the piston seal



CAUTION

Avoid touching the sealing surface in the pump head (17-4) with the tool.

Screw the special tool for the piston seal (31-1) with the narrow side just as far into the piston seal as the same can be removed.

5.5 High pressure pump

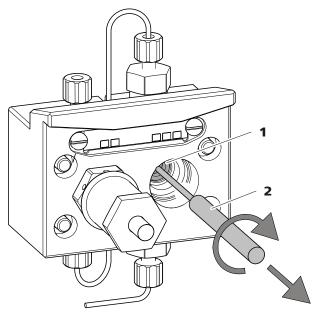


Figure 32 Removing the piston seal

1 Piston seal

Tool for piston seal Pin of the tool.

2 Inserting the new piston seal into the tool

Insert the new piston seal tightly by hand into the recess of the sleeve of the tool for the piston seal (31-2). The sealing springs must be visible from the outside.

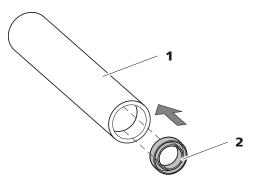


Figure 33 Inserting the piston seal into the tool

1 Tool for piston seal (6.2617.010)

Sleeve for inserting the new piston seal.

2 Piston seal

Order number: 6.2741.020

3 Inserting the new piston seal into the pump head

Guide the sleeve of the tool for the piston seal (31-2) with inserted piston seal into the pump head and press the seal with the wide end of the tool for the piston seal (31-1) into the pump head recess.

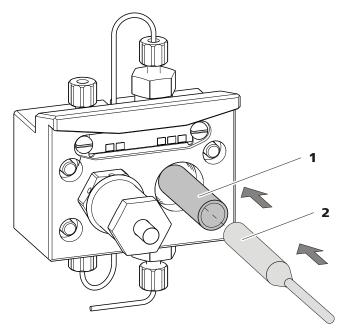


Figure 34 Inserting the piston seal into the pump head

4 Replacing the piston cartridge

Screw the assembled piston cartridge back into the pump head and tighten, first by hand, then additionally by approx. 15° with a wrench.

Cleaning the inlet valve and outlet valve

1 Removing valves

- Unscrew the connection capillary for the auxiliary piston (17-1) from the outlet valve holder.
- Unscrew the holders for the inlet and outlet valves and remove the valves (35-3) and (35-2).

5.5 High pressure pump

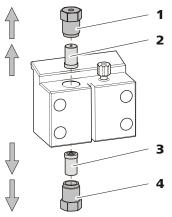


Figure 35 Removing valves

1 Outlet valve holder

Inlet valve

3

Order number: 6.2824.170

2 Outlet valve

Order number: 6.2824.160

4 Inlet valve holder

2 Cleaning undissected valve

Clean contaminated or blocked valves initially **without** dismantling them completely.

- Rinse the valve in eluent flow and counterflow direction using a spray bottle filled with ultrapure water, RBS solution or acetone.
- The rinsing effect is further increased through a short treatment (lasting for a maximum of 20 s) in an ultrasonic bath.



NOTICE

Longer lasting ultrasonic baths can damage the ruby ball of the valve.

Only if this cleaning is useless, dismantle the valves separately and clean the components.

3 Dismantling valve

Dismantle every valve separately.



NOTICE

For dismantling the valve the tool for valve cartridges (6.2617.020) is required.

- Place the valve with the seal faced downwards above the recess in the holder.
- Push the valve components out of the valve housing using the needle of the tool.

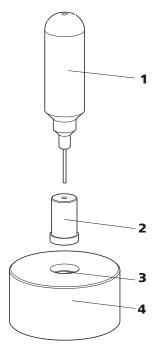


Figure 36 Dismantling valve

- Needle
 For pushing the valve components out of the valve housing.
- **3** Recess
 For collecting the valve components.

2 valve

4 Holder

The components of the valve are collected in the recess of the holder.



NOTICE

The components of the valve are very small. In order not to lose them, put the components into a dish.

• The inlet valve and the outlet valve consist of the same, just differently arranged components (see figure 37, page 68).

5.5 High pressure pump

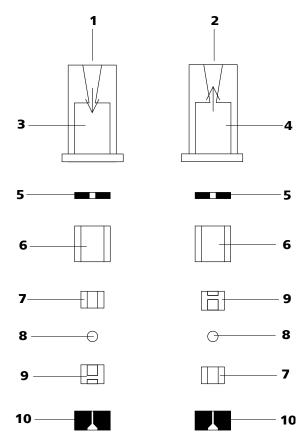


Figure 37 Components of the inlet valve and outlet valve

1	Inlet valve (6.2824.170)	2	Outlet valve (6.2824.160)
3	Inlet valve housing	4	Outlet valve housing
5	Sealing ring (black)	6	Sleeve
7	Sapphire sleeve The shiny side must point to ruby ball.	8	Ruby ball

4 Cleaning the components of the valve

Rinse the valve components with ultrapure water and/or acetone and dry with a lint-free cloth.

5 Reassembling the valve

Reassemble valve components according to figure 37, page 68.

- Insert the seal with the larger opening faced downwards into the recess of the tool.
- Lay the other valve components above another in the correct sequence (see figure 37, page 68).

5 Operation and maintenance

- Place over the valve housing and hold it tightly.
- By tilting the tool, the valve components slide into the valve housing.
- Press the seal by hand well on the valve housing.

6 Checking the flow direction

Rinse the valve in the direction of the arrow on the valve housing and check wether liquid is escaping on the other end.

If this is not the case, the valve has to be dismantled again and be reassembled correctly (see figure 37, page 68).

7 Inserting the valves back into the pump head



CAUTION

If by mistake, the inlet valve is mounted instead of the outlet valve, an extreme pressure builds up within the working cylinder, which can destroy the piston seal!

When inserting the valves, please take into account that the liquid is being pumped through the pump head from bottom to top.

- Insert the inlet valve into the inlet valve holder the way the seal is visible.
- Screw the inlet valve holder into the bottom of the pump head and tighten with a wrench (35-4).
- Insert the outlet valve into the outlet valve holder the way the seal is visible
- Screw the outlet valve holder into the top of the pump head and tighten with a wrench (35-1).

5.6 Inline filter -----

Mounting the pump head



NOTICE

To prevent the pump head from being positioned the wrong way, it is provided with different bore hole depths for the fastening bolts, i. e. a fastening bolt is longer than all others. The bore hole with the greatest depth must therefore be assigned to the longest bolt. If this is not the case, the pump will not function perfectly.


- 1 Mount the pump head on the pump again using the four fastening screws (17-**5**). Firmly tighten the screws with the hexagon key (6.2621.030).
- 2 Screw connection capillaries (17-1), (17-7) and (17-13) onto the pump head again.

5.6 **Inline filter**

5.6.1 **Maintenance**

The inline filter (6.2821.120) comprises the filter housing (38-2), the filter screw (38-4) and the filter (38-3). New filters (38-3) are available under the order number 6.2821.130 (10 items).

The filters (6.2821.130) (38-3) should be changed every 3 months (more frequently at higher backpressure).

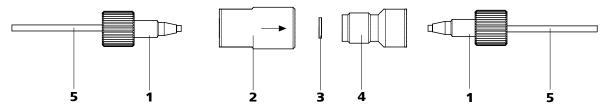

Change filters (of the inline filter) Figure 38

- PEEK pressure screws, short (6.2744.070)
- 3 Filter (6.2821.130) Packaging contains 10 items.
- 5 **Connection capillaries**

- 2 Filter housing Housing of the inline filter. Part of the 6.2821.120 accessories.
- Filter housing

Screw of the inline filter. Part of the 6.2821.120 accessories.

5 Operation and maintenance

Changing the filter

The flow must be stopped before changing the filter.

1 Removing the inline filter

• Unscrew the pressure screws (38-1) from the inline filter.

2 Unscrewing the filter screw

• Screw the filter screw (38-4) out of the filter housing (38-2) with the aid of two adjustable wrenches (6.2621.000).

3 Inserting the filter

- Remove the old filter (38-3) using tweezers.
- Place the new filter (38-3) flat in the filter housing (38-2) using tweezers.

4 Mounting filter screw

• Screw the filter screw (38-4) back into the filter housing (38-2) and tighten by hand. Then additionally tighten slightly with two adjustable wrenches (6.2621.000).

5 Remounting the inline filter

• Screw pressure screws (38-1) back onto the inline filter.

6 Rinsing the inline filter

- Dismantle the guard column (if present) and the separation column and replace with a coupling (6.2744.040).
- Rinse the instrument with eluent.

5.7 Inline sample preparation

To protect the separation column (see chapter 3.20, page 50) against foreign particles which can affect the separating efficiency, we recommend that all samples undergo a microfiltration (filter 0.45 μ m). The ultrafiltration cell can be used for **filtration** (see manual of the *IC Equipment for Ultrafiltration*).

Samples containing significant amounts of **gas** should be degassed. The sample degasser (see chapter 3.13, page 38) is used for degassing.

Matrix-loaded samples (e.g. blood, oil) should be prepared for the measurement by means of dialysis (see manual on the *IC Equipment for Dialysis*).

If the concentration of the sample is too high, the sample should be **diluted** before feeding (see documentation on the *IC Equipment for Sample Dilution*).

For the sample preparation methods **Neutralization** (replacement of e.g. Na⁺ with H⁺) and **cation exchange** (replacement of e.g. heavy metals with H⁺), a sample preparation module (SPM) is used.

For an overview of all Metrohm inline sample preparation methods go to the following website: http://misp.metrohm.com

5.8 Rinsing the sample path

Before a new sample can be measured, the sample path must be rinsed with it so that the measuring result is not falsified by the previous sample (**Sample carry-over**).

In the case of automated sample feeding, the rinsing time should be at least 3 times the **transfer time**.

The transfer time is the time required by the sample to flow from the sample vessel to the end of the sample loop. The transfer time depends on the pump capacity of the peristaltic pump or the Dosino used, the total capillary volume and the volume of the gas removed by the sample degasser (if used) - in other words the amount of gas in the sample.

5 Operation and maintenance

Ascertaining the transfer time

To ascertain the transfer time, proceed as follows:

1 Emptying the sample path

Pump air through the sample path (pump tubing, tubing connections, capillary in the degasser, sample loop) for several minutes until all liquid is displaced by the air.

2 Aspirating the sample and measuring time

Aspirate a sample typical for the later application and use a stop watch to measure the time required by the sample to travel from the sample vessel to the end of the sample loop.

The time measured corresponds to the "transfer time". The rinsing time should be at least 3 times the transfer time.

Checking the rinsing time

It is possible to determine whether the rinsing time is adequate via a direct measurement of the sample carry-over. Proceed as follows:

1 Preparing two samples

- **Sample A**: A typical sample for the application.
- Sample sample B: Ultrapure water.

2 Determining "Sample A"

Let "Sample A" pass through the sample path for the duration of the rinsing time, then inject and measure.

3 Determining "Sample B"

Let "Sample B" pass through the sample path for the duration of the rinsing time, then inject and measure.

4 Calculating the sample carry-over

The degree of the sample carry-over corresponds to the ratio of the peak areas of the measurement for sample B to the measurement for sample A. The lower the ratio, the lower the sample carry-over. This ratio can be modified by varying the rinsing time – thus allowing the rinsing time required for the application to be ascertained.

5.9 Sample degasser

5.9 Sample degasser

5.9.1 Operation

If you are working with sample degassing, the longer "transfer time" (see Ascertaining the transfer time, page 73) means that it is also necessary to rinse for longer (with the subsequent sample). The rinsing time should be at least 3 times the "transfer time" so as to minimize the carry-over effects. The "transfer time" depends on the pump capacity, the total capillary volume and the volume of the removed gas (i.e. the gas amount in the sample).



NOTICE

The rinsing time increases by at least 2 minutes when the sample degasser is used.

5.10 Injection valve

5.10.1 Protection

To prevent contamination of the injection valve, a 6.2821.120 inline filter (see chapter 3.11, page 36) should be mounted between the high pressure pump and the pulsation damper.

5.11 Conductivity detector

5.11.1 Maintenance



CAUTION

The conductivity detector must not be opened!



WARNING

When rinsing the detector, the pressure must not exceed **5 MPa**. In order to ensure this, the maximum pressure of the high pressure pump must be set to **5 MPa** in MagIC Net.

If the conductivity detector is blocked, you should first of all check whether the blockage results from the capillary ends being pressed together too firmly. In this case, shorten the detector input capillary (26-3) or the detector output capillary (27-2) by several millimeters.

If this does not help, the conductivity detector can be rinsed against the normal flow direction. For this, connect the high pressure pump to the detector output capillary (27-2) and rinse - the pressure may not exceed 5 MPa.

5.12 Separation column

5.12.1 Separating efficiency

Which analysis quality can be attained, depends to a great extent on the separating efficiency of the separation column used. The separating efficiency of the selected separation column must be sufficient for the analysis problems present. If difficulties occur, you should always first check the quality of the separation column by recording a standard chromatogram.

You can find detailed information on the separation columns available from Metrohm in the leaflet provided along with your separation column, in the **Metrohm IC-Column Program** (available via your Metrohm agent) or in the Internet at http://www.metrohm.com in the product area Ion chromatography. You can request free information on special IC applications in the corresponding "**Application Bulletins**" or "**Application Notes**", which are available in the Internet at http://www.metrohm.com in the Applications area or via the Metrohm agent responsible.

5.12.2 Protection

To protect the separation column against foreign particles, which can affect the separating efficiency, we recommend that both the eluent and the samples undergo a microfiltration (filter 0.45 μ m) before being aspirated via the aspiration filter (6.2821.090).

We recommend always to use a guard column (see chapter 3.19, page 49). This protects the separation column and considerably increases its service life. Information regarding which guard column is suitable for your separation column can be found in the **Metrohm IC Column Program** (which is available from your Metrohm agent), the leaflet provided along with your separation column, the product information on the separation column at http://www.metrohm.com (product area Ion Chromatography) or obtained directly from your agent.

The pulsation absorber (see chapter 3.12, page 37) must be installed in order to protect the column material from pressure concussion caused by injection.

5.12 Separation column

5.12.3 Storage

Always store the separation columns sealed and filled according to the data of the column manufacturer when not using them.

5.12.4 Regeneration



NOTICE

The regeneration is considered as the last measure, and not to be carried out regularly.

If the separating properties of the column have deteriorated, the column can be regenerated according to the specifications of the column manufacturer. In the case of separation columns available from Metrohm, the specification for regeneration can be found on the leaflet provided along with each column.

6 Troubleshooting

6 Troubleshooting

6.1 Problems and their solutions

Problem	Cause	Remedy	
Marked drop in pressure.	Leak in the system.	Check all capillary connections and seal leaks, if necessary (see chapter 3.5, page 15).	
The baseline has a large amount of noise.	High pressure pump – con- taminated pump valves.	Clean pump valves (see chapter 5.5.2, page 60).	
noise.	Eluent – Leakage in eluent path.	Check eluent path.	
	Eluent – Blockage in eluent path.	Check eluent path.	
	High pressure pump – defective piston seals.	Replace <i>(see chapter 5.5.2, page 60)</i> piston seals .	
	Pulsation damper not con- nected.	Connect the pulsation damper (see Chapter 3.12, Page 37).	
	Pulsation absorber not connected. or defective.	Connect pulsation absorber (see chapter 3.12, page 37). or replace it.	
The baseline is drifting.	Thermal equilibrium not yet attained.	Condition (see Chapter 4.2, Page 54) instrument with the column thermostat (see Chapter 3.15, Page 42) switched on .	
	Leak in the system.	Check all capillary connections and seal leaks, if necessary (see chapter 3.5, page 15).	
	Eluent – Evaporation of organic solvent in eluent.	 Check the eluent bottle cap (see figure 13, page 28). Stir the eluent. 	
The pressure in the system markedly increases.	Inline filter (6.2821.120) blocked.	Replace the filter (6.2821.130) (see chapter 5.6, page 70).	
mercases.	Conductivity detector blocked.	 Shorten capillary ends by several mm (see chapter 5.11.1, page 74). Rinse detector against the normal flow direction (see chapter 5.11.1, page 74). 	

Problem	Cause	Remedy		
	Guard column – blocked.	uard column – blocked. Replace guard column (see chapter 3.19, page 49).		
	Separation column – blocked.	 Regenerate separation column (see chapter 5.12.4, page 76). Replace separation column (see "Connecting and rinsing the separation column", page 51). Note: Samples should always be microfiltered 		
		(see chapter 5.7, page 72).		
	Injection valve – valve blocked.	Have the valve cleaned (by Metrohm service technicians).		
The retention times in the chromatogram have changed unexpectedly.	Separation column – diminished separating effi- ciency.	 Regenerate separation column (see chapter 5.12.4, page 76). Replace separation column (see "Connecting and rinsing the separation column", page 51). 		
	Eluent – Gas bubbles in eluent.	 Check the connectors of the eluent degasser (see chapter 3.9, page 30). Deaerate high-pressure pump (see chapter 3.10.2, page 34). 		
	High pressure pump – defective.	Request Metrohm Service.		
Peak areas are lower than expected.	Sample – leak in the sam- ple path.	Check the sample path.		
	Sample – blockage in the sample path.	Check the sample path.		
	Sample – sample loop not (completely) filled.	Prolong the sample transfer time.		
	Sample – gas bubbles in the sample.	Use sample degasser (see chapter 3.13, page 38).		
Data of the separa- tion column cannot	Column chip contami- nated.	Clean the contact surfaces of the column chip (with alcohol).		
be read.	Column chip defective.	 Save column configuration in MagIC Net™. Notify Metrohm Service. 		

6 Troubleshooting

Problem	Cause	Remedy
Individual peaks are greater than expected.	Sample – carry-over of the samples from previous measurement.	Rinse system longer between two samples.
The background conductivity is too high.	Wrong eluent.	Change the eluent (see chapter 5.4.2.3, page 59).
The retention times are poorly reprodu-	Eluent – Leakage in eluent path.	Check eluent path.
cible.	Eluent – Blockage in eluent path.	Check eluent path.
	Eluent – Gas bubbles in eluent.	 Check the connectors of the eluent degasser (see chapter 3.9, page 30). Deaerate high-pressure pump (see chapter 3.10.2, page 34).
Chromatograms have poor resolution	Separation column – diminished separating effi- ciency.	 Regenerate separation column (see chapter 5.12.4, page 76). Replace separation column (see "Connecting and rinsing the separation column", page 51).
Conductivity detector is not recognized in the software	No connection.	 Check the cable connection (27-1). Turn the instrument off and on again (after 15 seconds).
Extreme spread of Capillary connections the peaks in the chromatogram. Splitting (dual		Check connections (see chapter 3.5, page 15) (use PEEK capillaries with an internal diameter of 0.25 mm between the injection valve and detector).
peaks)	Guard column – diminished performance.	 Replace guard column (see chapter 3.19, page 49).
	Separation column – dead volume at column head.	 Install the separation column in reverse flow direction (if permitted by the leaflet) and rinse into a beaker. Replace separation column (see "Connecting and rinsing the separation column", page 51).
Precision problems - the measured values are highly scattered.	Sample – gas bubbles in the sample.	Use sample degasser (see chapter 3.13, page 38).

Problem	Cause	Remedy
	Injection valve – sample loop.	Check installation of the sample loop (see chapter 3.14.1, page 40).
	Sample – rinsing volume too low.	Increase rinsing time (see chapter 5.8, page 72).
	Injection valve – defective.	Request Metrohm Service.
Vacuum is not being built	Eluent degasser – Vacuum connector on the rear of the instrument not (tightly) sealed.	 Seal the Vacuum connector tightly with a threaded stopper (6.1446.040).

7 Technical specifications

7 Technical specifications

7.1 Reference conditions

The technical specifications listed in this chapter refers to the following

reference conditions:

Ambient tempera-

+25 °C (±3 °C)

ture

Instrument status > 40 minutes in operation (equilibrated)

7.2 Device

IC system ■ Metal-free IC system

• Compact system with modular design

• Up to two complete chromatographic systems in one housing

Material Painted polyurethane hard foam without CFCs, fire class V0

Operating pressure range

0...50 MPa (500 bar) high pressure pump0...35 MPa (350 bar) standard-PEEK system

Intelligent com-

iPump, iDetector, iColumn, MagIC Net

ponents

7.3 Leak sensor

Type Electronic, no calibration necessary

7.4 Ambient conditions

Operation

Ambient tem- +5...+45 °C

perature

Humidity 20...80 % relative humidity

Storage

Ambient tem- -20...+70 °C

perature

Transport

Ambient tem- -40...+70 °C

perature

7.5 Housing _____

Housing 7.5

Dimensions

Width 365 mm Height 642 mm Depth 380 mm

Material of base

Polyurethane hard foam (PUR) with flame retardation for fire class V0,

flask holder

tray, housing and CFC-free, painted

Operating ele-

ments

Indicators LED for power display

On/Off switch On the rear of the instrument

Eluent degasser 7.6

Material Fluoropolymer

Resistance to sol-

No restriction (except PFC)

vents

Time to establish

vacuum

< 60 s

High pressure pump 7.7

Туре Serial dual-piston pump

Intelligent pump head recognition

Chemically inert

Metal-free pump heads

Materials in contact with the eluent: PEEK, ZrO₂, PTFE/PE

Self-optimizing flow and pressure

Flow rate

Adjustable flow

0.001...20.0 mL/min

range

Flow increment 1 μL/min

Reproducibility

< 0.1 % deviation

of the eluent

flow

7 Technical specifications

Pressure range

Pump 0...50.0 MPa (0...500 bar)

< 1 %

Pump head 0...35.0 MPa (0...350 bar) (applies for the standard PEEK pump head)

Residual pulsa-

tion

Safety shutdown

Function Automatic shutdown upon reaching the pressure limit values

Maximum pressure limit Adjustable from 0.1...50 MPa (1...500 bar)

• The pump is automatically shut down at the first piston stroke above the maximum limit value

Minimum pressure limit

Adjustable from 0...49 MPa (0...490 bar)

• The shutdown mechanism is inactive at 0 MPa

The shutdown mechanism only becomes active 2 minutes after system start

• The pump is automatically shut down after 3 piston strokes below the minimum pressure limit

Gradient capacity Isocratic or gradient (extendable to quaternary)

Profile Step, linear, convex and concave Resolution < 1 nL/min flow increments

7.8 Sample degasser

Material Fluoropolymer

Resistance to sol-

vents

No restriction (except PFC)

Time to establish

vacuum

< 60 s

7.9 Injection valve

Actuator time typ.100 ms

Max. operating

pressure

Material

PEEK

35 MPa (350 bar)

7.10 Column thermostat

7.10 Column thermostat

Type Peltier technique thermostat for two intelligent separation columns

Adjustable tem-

0...+ 80 °C, in increments of 0.1 °C

perature range

Heating Ambient temperature +50 °C

Cooling Ambient temperature -20 °C

Temperature ± 0.2 °C

reproducibility

Stability < 0.05 °C

Heating up time < 30 minutes from 20 to 50 °C

Cooling time < 40 minutes from 50 to 20 °C

7.11 Conductivity measuring system

Type • Microprocessor-controlled Digital Signal Processing (DSP technol-

ogy)

Intelligent detector with 6 reference chromatograms

Measuring range 0...15000 μS/cm single range

Noise $< 0.1 \text{ nS at 1 } \mu\text{S/cm}$

Drift < 0.2 nS/cm per hour

Measuring rate 10 measurements per second for optimum results without filtering

Resolution 0.0047 nS/cm

Baseline Noise < 3 nS/cm typical without suppression

Conductivity detector

Cell volume 0.8 µL

Cell constant • Individual calibration data saved in the detector

adjustable range: 13.0...21.0 /cm

Electrodes Ring-shaped electrodes made from stainless steel

Materials in Chemically inert PCTFE

contact with

eluent

Maximum oper- 5.0 MPa (50 bar)

ating pressure

Cell tempera- 20...50 °C in increments of 5 °C

ture

7 Technical specifications

Temperature < 0.001 °C

stability

0...5 %/K adjustable, default 2.3 %/K

Temperature compensation

Heating up time < 30 minutes (40 °C)

7.12 Power connection

Required supply $100 - 240 \text{ V} \pm 10\%$ (autosensing)

voltage

quency

Required fre-

50 - 60 Hz \pm 3 Hz (autosensing)

Power consump-

■ 65 W for typical analysis application

tion

25 W standby (conductivity detector to 40 °C)

Power supply unit

• Up to 300 W maximum, electronically monitored

Internal fuse 3.15 A

7.13 Interfaces

USB

Entry 1 USB upstream, type B (for connection to the PC)

Exit 2 USB downstream, type A

MSB 2 MSB 8-pin Mini DIN (female) (for Dosino, stirrer, remote lines, etc.)



CAUTION

When connecting an instrument to the MSB connector you **must** switch off the 850 Professional IC.

Detector 2 15-pin high-density DSUB (female)

Column recogni-

tion

3 (including 2 in the column thermostat (see chapter 3.15, page 42))

Leak sensor 1 jack plug

Further connec-

.

■ 1 15-pin DSUB (female)

tions

7.14 Weight

7.14 Weight

1.850.3000 30.1 kg (without accessories)

1.850.9010 (con-

ductivity detector)

Transport carriage (rollers and han-

dle)

1.8 kg

2.3 kg (with accessories)

8 Accessories

8 Accessories

Up-to-date information on the scope of delivery and optional accessories for your product can be found on the Internet. You can download this information using the article number as follows:

Downloading the accessories list

- **1** Enter *https://www.metrohm.com/* into your Internet browser.
- 2 Enter the article number (e.g. **2.850.3000**) into the search field. The search result is displayed.
- Click on the product.

 Detailed information regarding the product is shown on various tabs.
- 4 On the **Included parts** tab, click on **Download the PDF**.

 The PDF file with the accessories data is created.



NOTICE

Once you have received your new product, we recommend downloading the accessories list from the Internet, printing it out and keeping it together with the manual for reference purposes.

Index

Index

(6.2821.130) Filter	70	Contamination		Filter (6.2821.130) 70	C
6.2821.090 aspiration filter	59	High pressure pump	59	Filter 6.2821.090	
		Valves in the high pressure		Aspiration filter 59	9
Α		pump	60	Flow fluctuations 60	C
Ambient conditions	81	Crystallization		Flow increment 82	2
Anion channel	. 7	High pressure pump	59	Flow range 82	2
Aspiration filter 6.2821.090	59	- ' ' '		Flow rate 82	2
Aspiration tubing for eluent	26	D		Frequency 85	5
		Deaerating			
В		High pressure pump	34	G	
Baseline		Purge valve		Gas 30, 38	3
Conditioning	55	Degasser		Guard column	
Unstable	60	Eluent degasser	30	Installation 49	9
Blockage		Sample degasser		Rinsing 50	C
Conductivity detector	74	Degassing		9	
Blood		Eluent	. 30	н	
		Detector		Handle 17	7
C		Cable connector	. 20	Heating	
Cable feed-throughs	23	Conductivity detector		see also "Column thermostat"	
Capillaries		Interface		42	2
Installation	15	Positioning		High pressure pump	
Capillary feed-throughs	23	Diagram		Installation 32	2
Carry-over		Dilution		Maintenance 59	
Cation channel	_	Dimensions		Protection 20, 59	9
Cleaning		Door		Technical specifications 82	
Valves of the high pressure		Drainage tubings	30	Tubing connection 32	
pump		Installation	21	Valves 68	
Column		mistanation	21	Housing 82	
see also "Separation column"	ı	E		Humidity 8	
		= Electrostatic charge	4		
Column recognition		Eluent	'	1	
Column thermostat		Aspirate	26	IC column	-
Column thermostat	84	Change		see also "Separation column"	
Installation		Production		50)
Computer connection		Eluent bottle	. 50	Initial installation 10	
Conditioning		Figure	29	Inject	•
Conductivity detector	33	Installation		Injection valve 42	2
*	45	Operation		Injection valve	-
Cell constant		Eluent degasser	. 33	Fill 42	2
Cell volume		Installation	20	Inject 42	
Maintenance				Installation 40, 83	
		Technical specifications		Maintenance	
Conductivity measuring system		Equilibration	. 54	Protection 74	
Technical data		Г		Inline filter 36	
Connect		Food throughs			
Power grid		Feed-throughs	22	Inline sample preparation 72	
To computer		Capillaries	. 23	Installation	
Connection		Fill	42	Conductivity detector	
Mains		Injection valve	42	Connections 45	
Connections		Filter	26	Connections	
Installation	15	see also "inline filter"	. 36	Drainage tubings 2°	1

Index

MSB 85 Pressure limit 83 Start-up 53 USB 85 Pressure range 83 Storage 81 Interfaces 85 Pressure range 83 Storage 81 Interface 85 Pressure range 83 Storage 81 Eurher connections 85 Pressure range 83 Storage 81 Euak sensor 85 Connection 15 Technical data Column thermostat 84 Leak sensor Pulsation absorber 60 Conductivity measuring system 84 Leak sallation 20 Purp head Detector 85 Technical specifications 81 Purp head Detector 85 Technical specifications 81 Purp head Detector 85 Technical specifications 81 Reference conditions	Eluent bottle 26	Overwiev of the instrument	Screws
High pressure pump	Eluent degasser 30	Front 6	Connection 15
Initial installation	Guard column 49		Separation column
Injection valve	High pressure pump 32	P	Installation 50
Leak sensor 20	Initial installation 10	Piston seal 60	Protection 37, 75
Pulsation absorber 37	Injection valve 40, 83	Pistons of the high pressure pump	Regeneration 76
Sample degasser 38	Leak sensor 20	60	Rinsing 52
Separation column	Pulsation absorber 37	Power connection 48, 85	Separating efficiency 75
Interface	Sample degasser 38	Power consumption 85	Storage 76
MSB 85 Pressure limit 83 Start-up 53 USB 85 Pressure range 83 Storage 81 Interfaces 85 Pressure screws Supply voltage 3,85 Further connections 85 Connection 15 Leak sensor 85 Protection Technical data Leak sensor 60 Pulsation 60 Conductivity measuring system Leak sensor 9 Installation 37 Technical data Conductivity measuring system Leak sensor 9 Installation 37 Technical specifications 84 Connection socket 9 Installation 37 Technical specifications 84 Interface 85 Maintenance 60 Eluent degaser 82 Technical specifications 81 Reference conditions 81 Reference conditions 81 Leak rightness 53.54 Reference conditions 81 Reference conditions 81 Reference conditions 81	Separation column 50	Power supply unit 85	Service 3, 56
USB	Interface	Precipitates 59	Shutting down 57
Interfaces	MSB 85	Pressure limit 83	Start-up 53
Further connections	USB 85	Pressure range 83	Storage 81
Leak sensor 85	Interfaces 85	Pressure screws	Supply voltage 3, 85
Injection valve	Further connections 85	Connection 15	
Leak	Leak sensor 85	Protection	Т
Leak Separation Separatio		Injection valve 74	Technical data
Leak	L	Inline filter 36	Column thermostat 84
Installation 37 Technical specifications 38 Detector 85 Maintenance 60 Eluent degasser 32 High pressure pump 82 Leak sensor 81 Reference conditions 82 Research 81 Reference conditions 81	Leak 60	Pulsation 60	Conductivity measuring system
Installation	Leak sensor	Pulsation absorber	84
Interface	Connection socket 9	Installation 37	Technical specifications
Technical specifications	Installation 20	Pump head	Detector 85
Leak-tightness 53, 54 Leaking piston seals 60 R Reference conditions 81 Reference conditions 82 Reference conditions 81 Reference conditions 82 Reference conditions 82 Reference conditions 83 Reference conditions 83 Reference conditions 84 Reference conditions 85 Reference conditions 85	Interface 85	Maintenance 60	Eluent degasser 82
Leaking piston seals 60 R Leak sensor 81 Loop Reference conditions 81 Reference conditions 81 see also "Sample loop" 42 Regeneration 56 Sample degasser 83 Mins connection Guard column 50 Separation column 50 see also "Column thermostat" Mains connection socket 8 Sample path 72 42 see also "Column thermostat" Mains connection socket 8 Sample path 72 42 see also "Column thermostat" Maintenance Separation column 52 Transfer time 73 Transfer time 73 Transport 81 81 81 81 81 81 81 82 82 82 83 83 84	Technical specifications 81	Purge valve32	High pressure pump 82
Leaking piston seals	Leak-tightness 53, 54	•	
Reference conditions Satisfactor Satis	Leaking piston seals 60	R	Leak sensor 81
Rinsing	Loop	Reference conditions 81	Reference conditions 81
Rinsing	see also "Sample loop" 42	Regeneration 56	Sample degasser 83
M Conductivity detector 74 Thermostat Mains connection Guard column 50 see also "Column thermostat" Mains connection socket 8 Sample path 72 42 Maintenance Separation column 52 Transfer time 73 Conductivity detector 74 Rinsing time 73 Transport 81 High pressure pump 59 Rise in pressure 59 Rollers 17 Injection valve 74 Rollers 17 Transport 81 Pump head 60 17 Transport locking screws 20 Material 82 5 Transport locking screws 20 Massuring range 84 Safety instructions 3 Installation 15 MPak Safety shutdown 83 Installation 15 MSB 85 Carry-over 72 USB 85 Connectors 9 Sample degasser V Vacuum pump Noise <td< td=""><td></td><td>Rinsing</td><td>· -</td></td<>		Rinsing	· -
Mains connection socket 8 Sample path 72 42 Maintenance Separation column 52 Transfer time 73 Conductivity detector 74 Rinsing time 73 Transport 81 High pressure pump 59 Rise in pressure 59 Rollers 17 Injection valve 74 Rollers 17 Transport locking screws 20 Pump head 60 Tubing 12 Tubing 12 Material 82 S Tubings 12 Measuring range 84 Safety instructions 3 Installation 15 MPak Safety shutdown 83 Installation 15 MSB Sample U USB 85 Connectors 9 Sample loop 42 Connectors 9 N Sample degasser V Vacuum pump Vacuum pump Protection 20 N Technical specifications 83 Valve Se	M	Conductivity detector 74	•
Maintenance Separation column 52 Transfer time 73 Conductivity detector 74 Rinsing time 73 Transport 81 High pressure pump 59 Rise in pressure 59 Rollers 17 Injection valve 74 Rollers 17 Transport locking screws 20 Pump head 60 Tubing 12 Material 82 Tubings 12 Measuring range 84 Safety instructions 3 Installation 15 MPak Safety shutdown 83 Installation 15 MSB Sample U USB 85 Connectors 9 Sample loop 42 Connectors 9 Noise 84 Installation 38 Vacuum pump Your under the present of the high pressure pump Oberation 72 Sample loop 42 see also "Injection valve" 40 Operation Sample path Valves of the high pressure pump 68	Mains connection	Guard column 50	see also "Column thermostat"
Maintenance Separation column 52 Transfer time 73 Conductivity detector 74 Rinsing time 73 Transport 81 High pressure pump 59 Rise in pressure 59 Rollers 17 Injection valve 74 Rollers 17 Transport locking screws 20 Pump head 60 Tubing 12 Material 82 Safety instructions 3 Installation 15 Measuring range 84 Safety shutdown 83 Installation 15 MPak Sample U USB 85 MSB 85 Carry-over 72 USB 85 Connectors 9 Sample loop 42 Connectors 9 Noise 84 Installation 38 Vacuum pump Operation 74 Protection 20 O Technical specifications 83 Valve Sample degasser 74 Sample loop	Mains connection socket 8	Sample path 72	42
High pressure pump 59	Maintenance		Transfer time 73
High pressure pump 59	Conductivity detector 74	Rinsing time 73	Transport 81
Injection valve			Rollers 17
Pump head 60 Material 82 Safety instructions 3 Tubings 12 Measuring range 84 Safety shutdown 83 MPak Safety shutdown 83 Holder 19 Sample U MSB 85 Carry-over 72 USB 85 Connectors 9 Sample loop 42 Connectors 9 Noise 84 Installation 38 Vacuum pump Noise 84 Installation 38 Vacuum pump Operation 74 Protection 20 Oil 72 Sample loop 42 see also "Injection valve" 40 Operation 38 Valves of the high pressure pump 68 Overview of the instrument 6 Sample preparation 72 68		Rollers 17	Transport locking screws 20
Material 82 S Tubings Measuring range 84 Safety instructions 3 MPak Safety shutdown 83 Holder 19 Sample MSB 85 Carry-over 72 Connectors 9 Sample loop 42 Connectors 9 N Sample degasser V Noise 84 Installation 38 Vacuum pump Operation 74 Protection 20 Oil 72 Sample loop 42 see also "Injection valve" 40 Operation 72 Sample path Valves of the high pressure pump 40 Operation of the instrument 74 Rinsing 72 72 68 Overview of the instrument 6 Sample preparation 72 72 68	Pump head 60		
Measuring range 84 Safety instructions 3 Installation 15 MPak Safety shutdown 83 Holder 19 Sample U MSB 85 Carry-over 72 USB 85 Connectors 9 Sample loop 42 Connectors 9 N Sample degasser V Noise 84 Installation 38 Vacuum pump Operation 74 Protection 20 Oil 72 Sample loop 42 see also "Injection valve" 40 Operation 5ample path Valves of the high pressure pump 68 Overview of the instrument 6 Sample preparation 72 68	Material 82	3	_
MPak Safety shutdown 83 Holder 19 Sample U MSB 85 Carry-over 72 USB 85 Connectors 9 Sample loop 42 Connectors 9 N Sample degasser V Noise 84 Installation 38 Vacuum pump Operation 74 Protection 20 O Technical specifications 83 Valve Operation 72 Sample loop 42 see also "Injection valve" 40 Operation Sample path Valves of the high pressure pump 68 Overview of the instrument 6 Sample preparation 72	Measuring range 84	Safety instructions 3	_
MSB 85 Carry-over 72 USB 85 Connectors 9 Sample loop 42 Connectors 9 N Sample degasser V Noise 84 Installation 38 Vacuum pump Operation 74 Protection 20 Valve Sample loop 42 see also "Injection valve" 40 Operation Sample path Valves of the high pressure pump 40 Sample degasser 74 Rinsing 72 72 68 Overview of the instrument 6 Sample preparation 72 72 68	MPak		
MSB 85 Carry-over 72 USB 85 Connectors 9 Sample loop 42 Connectors 9 Iransfer time 73 N Sample degasser V Noise 84 Installation 38 Vacuum pump Operation 74 Protection 20 Valve See also "Injection valve" 40 Operation Sample loop 42 see also "Injection valve" 40 Operation Sample path Valves of the high pressure pump 68 Overview of the instrument 6 Sample preparation 72	Holder 19	Sample	U
Connectors 9 Sample loop 42 Connectors 9 Inansfer time 73 73 73 73 73 74	MSB 85	Carry-over 72	USB 85
Transfer time 73 Noise 84 Installation 38 Vacuum pump Operation 74 Protection 20 O Technical specifications 83 Valve Oil 72 Sample loop 42 see also "Injection valve" 40 Operation Sample path Valves of the high pressure pump Sample degasser 74 Rinsing 72 68 Overview of the instrument 6 Sample preparation 72		Sample loop 42	Connectors 9
Noise 84 Installation 38 Vacuum pump Operation 74 Protection 20 O Technical specifications 83 Valve Oil 72 Sample loop 42 see also "Injection valve" 40 Operation Sample path Valves of the high pressure pump Sample degasser 74 Rinsing 72 Overview of the instrument 6 Sample preparation 72		Transfer time 73	
Operation 74 Protection 20 Oil 72 Sample loop 42 see also "Injection valve" 40 Operation Sample path Valves of the high pressure pump Sample degasser 74 Rinsing 72 Overview of the instrument 6 Sample preparation 72	N	Sample degasser	V
Operation 74 Protection 20 Oil 72 Sample loop 42 see also "Injection valve" 40 Operation Sample path Valves of the high pressure pump Sample degasser 74 Rinsing 72 Overview of the instrument 6 Sample preparation 72	Noise 84	Installation 38	Vacuum pump
OTechnical specifications83ValveOil			
Oil72Sample loop42see also "Injection valve"40OperationSample pathValves of the high pressure pumpSample degasser74Rinsing72Overview of the instrument6Sample preparation72	0	•	
OperationSample pathValves of the high pressure pumpSample degasser74Rinsing72Overview of the instrument6Sample preparation72	Oil 72	•	
Sample degasser	Operation		
Overview of the instrument 6 Sample preparation 72	'		
	Overview of the instrument 6	_	
	Rear 8	1 1 1 2 2 2 2 2 2 2 2 2 2 2 2 2 2 2 2 2	