917 Coulometer



Manual – Short Instructions 8.917.8002EN / v7 / 2024-10-30





info@metrohm.com www.metrohm.com

917 Coulometer

Manual - Short Instructions

Technical Communication Metrohm AG CH-9100 Herisau

This documentation is protected by copyright. All rights reserved.

This documentation is an original document.

This documentation has been prepared with great care. However, errors can never be entirely ruled out. Please send comments regarding possible errors to the address above.

Disclaimer

Deficiencies arising from circumstances that are not the responsibility of Metrohm, such as improper storage or improper use, etc., are expressly excluded from the warranty. Unauthorized modifications to the product (e.g. conversions or attachments) exclude any liability on the part of the manufacturer for resulting damage and its consequences. Instructions and notes in the Metrohm product documentation must be strictly followed. Otherwise, Metrohm's liability is excluded.

Melody for the BEEP command: excerpt from "En Altfrentsche", with kind permission of the Laseyer Quartett, Appenzell.

Table of contents

Table of contents

1	About thes	se short instructions	1
2	Introductio	n	2
	2.1 2.1.1 2.1.2 2.1.3	Instrument description Titration and measuring modes Connectors Intended use	2 3
	2.2 2.2.1	About the documentation	
3	Safety		5
	3.1	Intended use	5
	3.2	Responsibility of the operator	5
	3.3	Requirements for operating personnel	6
	3.4 3.4.1 3.4.2 3.4.3 3.4.4 3.4.5	Safety instructions General notes on safety Electrical safety Tubing and capillary connections Flammable solvents and chemicals Recycling and disposal	6 6 7
4	Overview o	of the instrument	9
	4.1	Front of the instrument	9
	4.2	Rear of the instrument	10
5	Installation	1	11
	5.1 5.1.1 5.1.2 5.1.3	Setting up the instrument Packaging Checks Location	11 11
	5.2 5.2.1 5.2.2 5.2.3	Setting up the titration cell for coulometry Inserting the titration cell Preparing the titration cell Mounting the addition and aspiration tube	12 12
	5.3	Connecting the coulometer to the power supply	17
	5.4 5.4.1 5.4.2 5.4.3	Connecting sensors Connecting generator electrodes Connecting indicator electrodes	18 20

Table of contents

	5.5	Connecting MSB devices	23
	5.5.1	Connecting a dosing device	24
	5.5.2	Connecting an additional stirrer or titration stand	25
	5.5.3	Connecting a Remote Box	25
	5.6	Connecting USB devices	26
	5.6.1	General	26
	5.6.2	Connecting a USB hub	26
	5.6.3	Connecting a printer	
	5.6.4	Connecting a balance	27
	5.6.5	Connecting a PC keyboard	28
	5.6.6	Connecting a barcode reader	
	5.6.7	Connecting the 885 Compact Oven SC	29
	5.7	Connecting the Coulometer to a network	30
6	Operation		31
	6.1	Switching the instrument on and off	31
	6.2	Fundamentals of operation	33
	6.2.1	Touch-sensitive screen	
	6.2.2	Display elements and controls	34
	6.2.3	Status display	35
	6.2.4	Entering text and numbers	36
7	Maintenan	ce	38

Table of figures

Table of figures

Figure 1	Front	9
Figure 2	Rear	10
Figure 3	Inserting the titration cell	12
Figure 4	Filling the adsorber tube	13
Figure 5	Equipping the titration cell	
Figure 6	Mounting the addition and aspiration tube	16
Figure 7	Connecting the power supply unit	17
Figure 8	Unscrewing the protective cap from the generator electrode	19
Figure 9	Screwing the electrode cable to the generator electrode	19
Figure 10	Connecting a generator electrode	20
Figure 11	Unscrewing the protective cap from the indicator electrode	20
Figure 12	Screwing the electrode cable to the indicator electrode	21
Figure 13	Connecting an indicator electrode	21
Figure 14	MSB connections	
Figure 15	MSB connector	24

1 About these short instructions

1 About these short instructions

This short instruction manual contains important information concerning the 917 Coulometer. In addition to an introduction, safety instructions and an overview of the instrument, you will also find information about the installation of the coulometer.

Additional information concerning the instrument can be found at http://products.metrohm.com.

2.1 Instrument description

2 Introduction

2.1 Instrument description

The 917 Coulometer is a compact titration system for coulometric Karl Fischer titration. This titrator combines in a single device the touch-sensitive color monitor for convenient and efficient operation, the magnetic stirrer and the membrane pump for adding and aspirating solvents. The upper side of the housing offers space for the titration cell. Thanks to its compact construction, you can use the 917 Coulometer in a small space as a stand-alone titrator.

You manage reagents, sensors, methods, etc. conveniently in the internal memory of the coulometer. You can also save your files externally through the USB connector, e.g. on a USB flash drive. On this storage medium you can not only store your methods and determinations, but also create a backup together with all of the data and settings of your system.

The integrated Ethernet connector is available to you should you wish to connect your 917 Coulometer to a network. The network connection offers you the following advantages:

- Saving data to a computer within the network
- Printing reports on a network printer
- Sending displayed messages as e-mails

2.1.1 Titration and measuring modes

The 917 Coulometer supports the following titration and measuring modes.

KFC

Coulometric water content determination according to Karl Fischer with voltametric measurement.

BRC

Coulometric bromine index determination. Determining the amount of double bonds in e.g. mineral oils.

MEAS

Measuring modes:

- Ipol (voltametric measurement with selectable polarization current)
- Upol (amperometric measurement with selectable polarization voltage)
- T (temperature measurement)

2 •••••• 917 Coulometer

2 Introduction

2.1.2 Connectors

The 917 Coulometer is equipped with the following connectors:

Power connection

For connecting to the power grid using the power supply unit provided

Two MSB connectors (Metrohm Serial Bus)

For connecting dosing devices, one additional stirrer or a Remote Box.

USB connector

For connecting peripheral devices (printer, PC keyboard, etc.), a USB flash drive, a Sample Processor or a USB hub.

Sensor connectors

One connector each for:

- Generator electrode
- Indicator electrode
- Temperature sensor (Pt1000 or NTC)

Ethernet connector

For connecting the Coulometer to a network.

Two connection nipples

For connecting tubing for replacing reagent using the built-in pump.

2.1.3 Intended use

The 917 Coulometer is designed for usage as a titrator in analytical laboratories. Its main application field is coulometric Karl Fischer titration.

This instrument is suitable for processing chemicals and flammable samples. Therefore, the use of the instrument requires the user to have basic knowledge and experience in handling toxic and caustic substances. Knowledge regarding the application of fire prevention measures prescribed for laboratories is also mandatory.

2.2 About the documentation



CAUTION

Please read through this documentation carefully before putting the instrument into operation. The documentation contains information and warnings which the user must follow in order to ensure safe operation of the instrument.

2.2 About the documentation

2.2.1 Symbols and conventions

The following symbols and formatting may appear in this documentation:

(5- 12)	Cross-reference to figure legend
	The first number refers to the figure number, the second to the instrument part in the figure.
1	Instruction step
	Perform the steps one after the other.
Method	Dialog text, parameter in the software
File ► New	Menu or menu item
[Continue]	Button or key
$\overline{\wedge}$	WARNING
	This symbol draws attention to a possible life-threat- ening hazard or risk of injury.
	WARNING
7	This symbol draws attention to a possible hazard due to electrical current.
	WARNING
	This symbol draws attention to a possible hazard due to heat or hot instrument parts.
	WARNING
	This symbol draws attention to a possible biological hazard.
	WARNING
深	Warning of optical radiation
A	CAUTION
	This symbol draws attention to possible damage to instruments or instrument parts.
i	NOTICE
	This symbol highlights additional information and tips.

4 ••••• 917 Coulometer

3 Safety

3 Safety

3.1 Intended use

The 917 Coulometer is designed for usage as a titrator in analytical laboratories. Its main area of use is coulometric Karl Fischer titration.

This device is suitable for processing chemicals and flammable samples. Therefore, the use of the device requires the user to have basic knowledge and experience in handling toxic and caustic substances. Knowledge of the application of the fire prevention measures prescribed for laboratories is also mandatory.

3.2 Responsibility of the operator

The operator must ensure that basic regulations on occupational safety and accident prevention in chemical laboratories are observed. The operator has the following responsibilities:

- Instruct personnel in the safe handling of the product.
- Train personnel in the use of the product according to the user documentation (e.g. install, operate, clean, eliminate faults).
- Train staff on basic occupational safety and accident prevention regulations.
- Provide personal protective equipment (e.g. protective glasses, gloves).
- Provide suitable tools and equipment to carry out the work safely.

The product may be used only when it is in perfect condition. The following measures are required to ensure the safe operation of the product:

- Check the condition of the product before use.
- Remedy defects and malfunctions immediately.
- Maintain and clean the product regularly.

3.3 Requirements for operating personnel

Only qualified personnel may operate the product. Qualified personnel are persons who meet the following requirements:

- Basic regulations on occupational safety and accident prevention for chemical laboratories are known and complied with.
- Knowledge of handling hazardous chemicals is present. Personnel have the ability to recognize and avoid potential dangers.
- Knowledge regarding the application of fire prevention measures for laboratories is available.
- Safety-relevant information is communicated and understood. The personnel can operate the product safely.
- The user documentation has been read and understood. The personnel operate the product according to the instructions in the user documentation.

3.4 Safety instructions

3.4.1 General notes on safety



WARNING

Operate this instrument only according to the information contained in this documentation.

This instrument left the factory in a flawless state in terms of technical safety. To maintain this state and ensure non-hazardous operation of the instrument, the following instructions must be observed carefully.

3.4.2 Electrical safety

The electrical safety when working with the instrument is ensured as part of the international standard IEC 61010.



WARNING

Only personnel qualified by Metrohm are authorized to carry out service work on electronic components.

6 •••••• 917 Coulometer

3 Safety



WARNING

Never open the housing of the instrument. The instrument could be damaged by this. There is also a risk of serious injury if live components are touched.

There are no parts inside the housing which can be serviced or replaced by the user.

Supply voltage



WARNING

An incorrect supply voltage can damage the instrument.

Only operate this instrument with a supply voltage specified for it (see rear panel of the instrument).

Protection against electrostatic charges



WARNING

Electronic components are sensitive to electrostatic charges and can be destroyed by discharges.

Do not fail to pull the power cord out of the power socket before you set up or disconnect electrical plug connections at the rear of the instrument.

3.4.3 Tubing and capillary connections



CAUTION

Leaks in tubing and capillary connections are a safety risk. Tighten all connections well by hand. Avoid applying excessive force to tubing connections. Damaged tubing ends lead to leakage. Appropriate tools can be used to loosen connections.

Check the connections regularly for leakage. If the instrument is used mainly in unattended operation, then weekly inspections are mandatory.

3.4 Safety instructions

3.4.4 Flammable solvents and chemicals



WARNING

All relevant safety measures are to be observed when working with flammable solvents and chemicals.

- Set up the instrument in a well-ventilated location (e.g. fume cupboard).
- Keep all sources of flame far from the workplace.
- Clean up spilled liquids and solids immediately.
- Follow the safety instructions of the chemical manufacturer.

3.4.5 Recycling and disposal



Properly dispose of chemicals and of the product to reduce negative effects on the environment and public health. Local authorities, waste disposal companies or dealers provide more detailed information on disposal. Observe the WEEE EU directive (WEEE = Waste Electrical and Electronic Equipment) for the proper disposal of waste electronic equipment within the European Union.

8 ------ 917 Coulometer

4 Overview of the instrument

4 Overview of the instrument

4.1 Front of the instrument

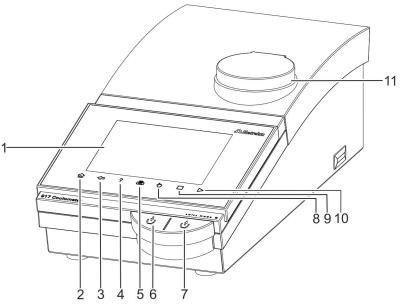


Figure 1 Front

1 Display

Touch screen.

3 Fixed key [Back]

Saves the entry and opens the next-higher dialog page.

5 Fixed key [Print]

Opens the print dialog.

7 Key

Pressing the key aspirates air out of the aspiration bottle. The vacuum in the aspiration bottle aspirates the liquid out of the titration cell and into the aspiration bottle.

9 Fixed key [STOP]

Cancels the running determination.

11 Magnetic stirrer

For placing the titration vessel holder.

2 Fixed key [Home]

Opens the main dialog.

4 Fixed key [Help]

Opens the online help for the dialog displayed.

6 Key

Pressing the key pumps air into the solvent bottle. The overpressure in the solvent bottle pushes solvent into the titration cell.

8 Fixed key [Manual]

Opens the manual control.

10 Fixed key [START]

Starts a determination.

4.2 Rear of the instrument

4.2 Rear of the instrument

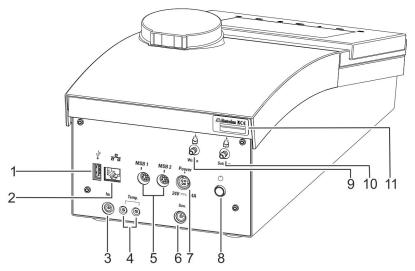


Figure 2 Rear

1 USB connector (type A)

For connecting printers, USB flash drives, USB hubs, 885 Compact Oven SC, etc.

- 3 Indicator electrode connector (Ind.)
 For connecting the indicator electrode.
- MSB connector (MSB 1 and MSB 2)
 Metrohm Serial Bus. For connecting external dosing devices, stirrers or a Remote Box.
 Mini DIN, 8-pin.
- **7** Power socket (Power)

For connecting the external power supply unit.

- **9 Connection nipple for PVC tubing**For aspirating the contents of the titration cell.
- 11 Type plate

Contains the serial number.

- Ethernet connector (RJ-45)
 For connecting to a network.
- **4** Temperature sensor connector (Temp.)
 For connecting temperature sensors (Pt1000 or NTC). Two B sockets, 2 mm.
- **Generator electrode connector (Gen.)** For connecting the generator electrode.
- 8 Power switch

Switches the instrument on/off.

10 Connection nipple for PVC tubing For aspirating solvent.

10 ----- 917 Coulometer

5 Installation

5.1 Setting up the instrument

5.1.1 Packaging

The instrument is supplied in protective packaging together with the separately packed accessories. Keep this packaging, as only this ensures safe transportation of the instrument.

5.1.2 Checks

Immediately after receipt, check whether the shipment has arrived complete and without damage by comparing it with the delivery note.

5.1.3 Location

The instrument has been developed for operation indoors and may not be used in explosive environments.

Place the instrument in a location of the laboratory which is suitable for operation and free of vibrations and which provides protection against corrosive atmosphere and contamination by chemicals.

The instrument should be protected against excessive temperature fluctuations and direct sunlight.

5.2 Setting up the titration cell for coulometry

5.2.1 Inserting the titration cell

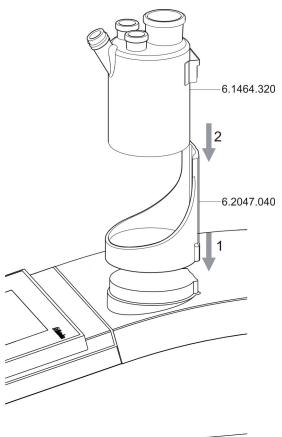


Figure 3 Inserting the titration cell

Set up the titration cell on the magnetic stirrer as follows:

1 Attach the 6.2047.040 titration vessel holder to the magnetic stirrer.

2 Insert the 6.1464.320 titration cell into the titration vessel holder.

5.2.2 Preparing the titration cell

Filling the adsorber tube

Before setting up the titration cell, the 6.1403.030 adsorber tube has to be filled with 6.2811.000 molecular sieve. Proceed as follows:

12 ••••• 917 Coulometer

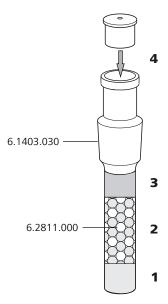


Figure 4 Filling the adsorber tube

- 1 Insert a small cotton plug into the bottom of the adsorber tube. Do not pack the cotton too tightly.
- **2** Fill the molecular sieve up to ¾ of the height.
- **3** Place a small cotton plug on the molecular sieve. Do not pack the cotton too tightly.
- **4** Seal the adsorber tube with the appropriate cover.



NOTE

Note that the molecular sieve must be replaced at regular intervals. Each time you refill the adsorber tube with molecular sieve, you can, for example, write the date directly on the adsorber tube.

6.0341.100 5 4 6.0345.100 6.0344.100 6.2713.000 6.2701.040 6.2701.040 6.1448.020 6.1448.020 6.1464.320 6.1903.030

Equipping the titration cell

Figure 5 Equipping the titration cell

Equip the titration cell as follows:

- 1 Place the 6.1903.030 stirring bar in the titration cell.
- 2 Cut the 6.2713.0x0 ground-joint sleeves to the correct length and attach them to the ground joints of the inserts (electrodes, adsorber tube, etc.).
 - Take care to ensure that the edges of the ground-joint sleeves are cut to size cleanly and that there are no fringes. The ground-joint sleeves must not protrude at the lower edge of the ground-joint opening.
- **3** Insert the 6.1403.030 adsorber tube into the generator electrode.

14 •••••• 917 Coulometer

- 4 Insert the 6.0345.100 generator electrode without diaphragm or the 6.0344.100 generator electrode with diaphragm together with the adsorber tube into the large ground-joint opening at the rear.
- Insert the 6.0341.100 indicator electrode into the left ground-joint opening.
- Place the 6.1448.020 septum on the front opening of the titration cell and screw it shut with the 6.2701.040 screw cap.

 Tighten the screw cap only enough so that it seals. The septum must not bend.

Filling the titration cell (generator electrode without diaphragm)

Proceed as follows when using a generator electrode without a diaphragm:

- **1** Fill approximately 100 mL of reagent into the titration cell using the 6.2738.000 funnel.
- 2 Close the remaining ground-joint opening on the right with the 6.1437.000 ground-joint stopper (with ground-joint sleeve attached).

Filling the titration cell (generator electrode with diaphragm)

Proceed as follows when using a generator electrode with a diaphragm:

- 1 Fill approximately 5 mL of catholyte into the generator electrode.
- Fill approximately 100 mL of anolyte into the titration cell using the 6.2738.000 funnel. The level of the anolyte should be roughly 1 2 mm above the level of the catholyte.
- Close the remaining ground-joint opening on the right with the 6.1437.000 ground-joint stopper (with ground-joint sleeve attached).

5.2.3 Mounting the addition and aspiration tube

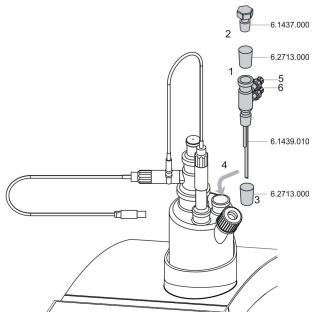


Figure 6 Mounting the addition and aspiration tube

Insert the addition and aspiration tube as follows into the titration cell:

- Attach the 6.2713.000 ground-joint sleeve that has been cut to size to the ground joint of the 6.1437.000 stopper.
- 2 Insert the stopper into the 6.1439.010 addition and aspiration tube.
- Attach the 6.2713.000 ground-joint sleeve that has been cut to size to the ground joint of the addition and aspiration tube.
- 4 Insert this assembly into the ground-joint opening.
- **5** Connect the tubing for the reagent addition at the upper connector of the addition and aspiration tube (5).
- 6 Connect the tubing for the aspiration of the titration cell at the lower connector of the addition and aspiration tube (6).

Details regarding how to connect the addition tubing and the aspiration tubing can be found in the tutorial for the 917 Coulometer.

16 ----- 917 Coulometer

5.3 Connecting the coulometer to the power supply

Connecting the power supply unit

The 917 Coulometer has an external power supply unit for a 24 V power supply (DC). This is connected to the power socket of the coulometer.



WARNING

An incorrect supply voltage can damage the instrument.

Operate the instrument only with the supply voltage specified for it. Use the supplied power supply unit exclusively, see chapter *Technical specifications*.

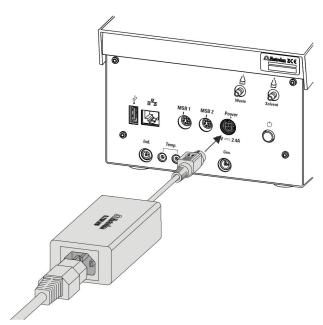


Figure 7 Connecting the power supply unit

Proceed as follows:

Connect the plug of the external power supply unit with the power socket of the coulometer.

5.4 Connecting sensors



NOTE

The plug of the power supply unit is protected against accidental disconnection of the cable by means of a pull-out protection feature. If you wish to pull out the plug, you first need to pull back the outer plug sleeve marked with arrows.

2 Connect the power cord with the external power supply unit of the coulometer and with the power grid.



CAUTION

Switch off the Coulometer correctly by pressing the power switch before you disconnect the electricity supply. If this is not done, then there is a danger of data loss.

5.4 Connecting sensors

The measuring interface contains the following measuring inputs:

- **Gen.** for a generator electrode
- Ind. for a double Pt electrode
- Temp. for a temperature sensor of the Pt1000 or NTC type



CAUTION

Under all circumstances, avoid mixing up the electrode cable from the indicator electrode with the one from the generator electrode. Mark the screw heads of the cables accordingly.

5.4.1 Connecting generator electrodes

Screwing the electrode cable to the generator electrode

1 Unscrew the protective cap from the generator electrode.

18 ----- 917 Coulometer

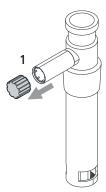


Figure 8 Unscrewing the protective cap from the generator electrode

2 Tighten the 6.2104.120 electrode cable to the generator electrode.

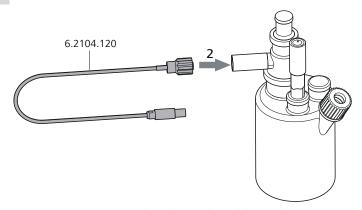


Figure 9 Screwing the electrode cable to the generator electrode

Connecting the electrode cable to the coulometer

1 Plug the electrode plug into the **Gen.** socket of the coulometer.

5.4 Connecting sensors

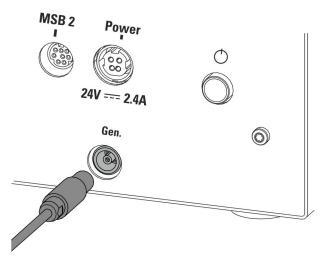


Figure 10 Connecting a generator electrode



NOTE

The electrode cable is protected against accidental disconnection of the cable by means of a pull-out protection. If you wish to pull out the plug again, you first need to pull back the outer plug sleeve.

5.4.2 Connecting indicator electrodes

Screwing the electrode cable to the indicator electrode

1 Unscrew the protective cap from the indicator electrode.

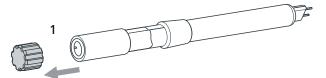


Figure 11 Unscrewing the protective cap from the indicator electrode

2 Tighten the 6.2104.020 electrode cable to the indicator electrode.

20 -----917 Coulometer

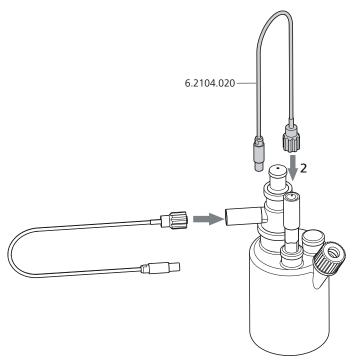


Figure 12 Screwing the electrode cable to the indicator electrode

Connecting the electrode cable to the coulometer

1 Plug the electrode plug into the **Ind.** socket of the coulometer.

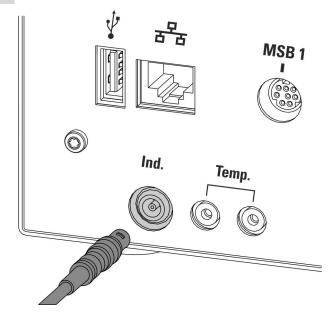


Figure 13 Connecting an indicator electrode

5.4 Connecting sensors



NOTE


The electrode cable is protected against accidental disconnection of the cable by means of a pull-out protection. If you wish to pull out the plug again, you first need to pull back the outer plug sleeve.

5.4.3 Connecting a temperature sensor

A temperature sensor of the Pt1000 or NTC type can be connected to the **Temp.** connector.

Connect the temperature sensor as follows:

1 Insert the plugs of the temperature sensor into the **Temp.** sockets of the coulometer.

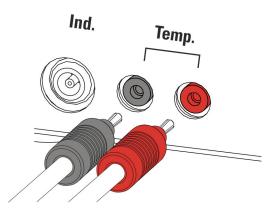



NOTE

Always insert the red plug into the red socket. This is the only way that shielding against electrical interference can be ensured.

22 •••••• 917 Coulometer

5.5 Connecting MSB devices

In order to connect MSB devices, e.g. dosing device or Remote Box, the Coulometer has two connectors at what is referred to as the *Metrohm Serial Bus* (MSB). Various peripheral devices can be connected in sequence (daisy chain) at a single MSB connector (8-pin Mini DIN socket) and be controlled simultaneously by the Coulometer. In addition to the connection cable, stirrers and the Remote Box are each equipped with their own MSB socket for this purpose.



NOTE

The additional stirrer must not be connected to the MSB 1!

The following figure provides an overview of the instruments that can be connected to an MSB socket, along with a number of different cabling variations.

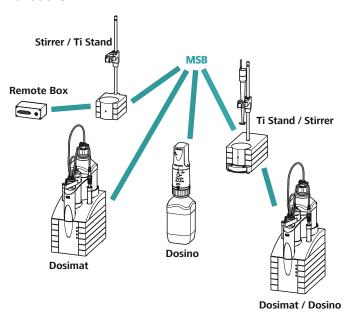


Figure 14 MSB connections



NOTE

When connecting MSB devices together, the following must be observed:

- No additional stirrer may be connected to the MSB 1!
- Only one device of the same type may be used at a single MSB connector at one time.

• When making the connection, take care to ensure that the flat part of the MSB plug marked with arrows is pointing in the direction of the marking on the MSB connector (see figure 15, page 24).

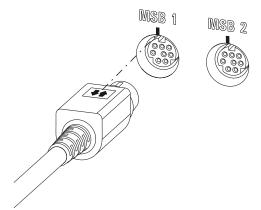


Figure 15 MSB connector



CAUTION

Switch off the Coulometer before you plug in MSB devices. When it is switched on, the Coulometer automatically recognizes which device is connected to which MSB connector. The connected MSB devices are entered automatically in the device manager.

MSB connections can be extended with the 6.2151.010 cable. The maximum connection length permitted is 6 m.

5.5.1 Connecting a dosing device

You can connect two dosing devices to the Coulometer.

The types of dosing devices that are supported are:

- 800 Dosino
- 805 Dosimat

Proceed as follows:

1 Switch off the Coulometer.

24 •••••• 917 Coulometer

2 Connect the dosing device connection cable to an MSB connector on the rear side of the Coulometer.

3 Switch on the Coulometer.

5.5.2 Connecting an additional stirrer or titration stand

Proceed as follows:

- **1** Switch off the Coulometer.
- **2** Connect the connection cable of the magnetic stirrer or of the titration stand to MSB 2 on the rear of the Coulometer.
- **3** Switch on the Coulometer.

5.5.3 Connecting a Remote Box

Instruments that are controlled via remote lines and/or that send control signals via remote lines can be connected via the 6.2148.010 Remote Box. In addition to Metrohm, other instrument manufacturers also use similar connectors that make it possible to connect different instruments together. These interfaces are also frequently given the designations "TTL Logic", "I/O Control" or "Relay Control" and they generally have a signal level of 5 volts.

Control signals are understood to be electrical line statuses or brief (> 200 ms) electrical pulses which display the operational state of an instrument or which trigger or report an event. Sequences on a variety of instruments can thus be coordinated in a single complex automation system. However, no exchange of data is possible.

Proceed as follows:

- 1 Switch off the Coulometer.
- **2** Connect the Remote Box connection cable to an MSB connector on the rear side of the Coulometer.
- **3** Switch on the Coulometer.

You will find precise information concerning the pin assignment of the interface on the Remote Box in the Appendix of the more detailed manual.

5.6 Connecting USB devices

5.6 Connecting USB devices

5.6.1 General

The 917 Coulometer has a USB connector (type A socket) for peripheral devices with a USB interface and for the 885 Compact Oven SC. If you wish to connect more than one device to the USB, you can use a commercially available USB hub.



NOTE

We recommend that the Coulometer be switched off while you set up or disconnect connections between the devices.

5.6.2 Connecting a USB hub

Use a USB hub with its own power supply.

Connect the USB hub as follows:

With the help of the 6.2151.030 cable (length 0.6 m) or the 6.2151.020 cable (length 1.8 m), connect the USB connector of the Coulometer (Type A) with the USB connector of the hub (Type B, see manual for the USB hub).

The USB hub is recognized automatically.

5.6.3 Connecting a printer

Printers connected to the 917 Coulometer must meet the following requirements:

- Printer languages: HP-PCL, HP-PCL-GUI, Canon BJL Commands or Epson ESC P/2
- Paper size: A4 or Letter.

Connect the printer as follows:

- With the aid of the 6.2151.020 cable, connect the USB connector of the Coulometer (type A) with the USB connector of the printer (type B, see manual for the printer).
- **2** Configure the printer in the device manager of the Coulometer.

26 ------

5.6.4 Connecting a balance

If you want to connect a balance to the Coulometer you need a USB/RS-232 adapter (6.2148.050).

The following table offers an overview of the balances that you can use together with the Coulometer and of which cable you will need for connection to the RS-232 interface:

Balance	Cables
AND ER, FR, FX with RS-232 interface (OP-03)	6.2125.020 + 6.2125.010
Mettler AB, AG, PR (LC-RS9)	In the scope of delivery for the balance
Mettler AM, PM, PE with interface	6.2146.020 + 6.2125.010
option 016 or Mettler AJ, PJ with interface	Also from Mettler: ME 47473 adapter and either ME 42500 hand switch or ME 46278 foot
option 018	switch
Mettler AT	6.2146.020 + 6.2125.010
	Also from Mettler: ME 42500 hand switch or ME 46278 foot switch
Mettler AX, MX, UMX, PG, AB-S, PB-S, XP, XS	6.2134.120
Mettler AE with interface option	6.2125.020 + 6.2125.010
011 or 012	Also from Mettler: ME 42500 hand switch or ME 46278 foot switch
Ohaus Voyager, Explorer, Analytical Plus	Cable AS017-09 from Ohaus
Precisa balances with RS-232-C interface	6.2125.080 + 6.2125.010
Sartorius MP8, MC, LA, Genius, Cubis	6.2134.060
Shimadzu BX, BW	6.2125.080 + 6.2125.010

Connect the balance as follows:

1 Connect the USB plug of the USB/RS-232 adapter with the USB connector of the Coulometer (Type A).

- The USB/RS-232 adapter will be recognized automatically and entered in the device manager of the Coulometer.
- **2** Connect the RS-232 interface of the USB/RS-232 adapter with the RS-232 interface of the balance (see table for cable).
- **3** Switch on the balance.
- **4** If necessary, switch on the RS-232 interface of the balance.
- **5** Configure the RS-232 interface of the USB/RS-232 adapter in the device manager of the Coulometer .
- **6** Enter and configure the balance in the device manager of the Coulometer.
- **7** Make sure that the parameters of the USB/RS-232 adapter configured in the device manager match those of the balance.

5.6.5 Connecting a PC keyboard

The PC keyboard is used as an aid for text and numerical input.

Connect the PC keyboard as follows:

- 1 Connect the USB plug of the keyboard with the USB connector of the Coulometer (Type A).
- **2** Enter and configure the keyboard in the device manager of the Coulometer.

5.6.6 Connecting a barcode reader

The barcode reader is used as an aid for text and numerical input. You can connect a barcode reader with USB interface.

Connect the barcode reader as follows:

1 Connect the USB plug of the barcode reader with the USB connector of the Coulometer (Type A).

28 •••••• 917 Coulometer

2 Enter and configure the barcode reader in the device manager.

Settings on the barcode reader:

Program the barcode reader as follows (also see manual for the barcode reader):

- **1** Switch the barcode reader to programming mode.
- **2** Specify the desired layout for the keyboard (USA, Germany, France, Spain, German-speaking Switzerland).

This setting must match the setting in the device manager.

- Make sure that the barcode reader is set in such a way that Ctrl characters (ASCII 00 to 31) can be sent.
- Program the barcode reader in such a way that the ASCII character 02 (STX or Ctrl B) is sent as the first character. This first character is normally referred to as the "Preamble" or "Prefix Code".
- Program the barcode reader in such a way that the ASCII character 04 (EOT or Ctrl D) is sent as the last character. This last character is normally referred to as the "Postamble", "Record Suffix" or "Postfix Code".
- **6** Exit the programming mode.

5.6.7 Connecting the 885 Compact Oven SC



NOTE

In order for the 885 Compact Oven SC to be controlled with the 917 Coulometer, the following firmware versions must be installed on the instruments as a minimum requirement:

- 917 Coulometer 5.917.0025
- 885 Compact Oven SC 5.885.0013

The 885 Compact Oven SC is connected via the USB interface. The 6.2151.110 USB cable is required for this.

If the USB interface on the 917 Coulometer is also needed for other USB devices (USB flash drive, USB keyboard, USB/RS-232 adapter for balance), use a USB hub with its own power supply.

Connect the 885 Compact Oven SC as follows:

- **1** Switch off the 917 Coulometer.
- **2** Connect the 885 Compact Oven SC to the Coulometer using a USB cable.

- **3** Switch on the 885 Compact Oven SC.
- **4** Switch on the 917 Coulometer.



NOTE

When the 917 Coulometer is started, the 885 Compact Oven SC is added to the device manager and the settings from the device manager are transferred to the instrument.

The switch-on sequence must be strictly observed with the 885 Compact Oven SC.

5 Configure the 885 Compact Oven SC in the device manager.

5.7 Connecting the Coulometer to a network

The 917 Coulometer has a network connection (Ethernet). This can be used to integrate your Coulometer in your network. You can, for example, store data on a PC within the network or print reports on a network printer. You will find information as to which settings are necessary for the network connection in the *Device manager* chapter contained in the more detailed manual.

Connecting the Coulometer to a network

30 ••••••

6 Operation

6 Operation

6.1 Switching the instrument on and off

Switching on the instrument



CAUTION

Peripheral devices (e.g., printers, 885 Compact Oven SC, etc) must be connected and switched on before you switch on the 917 Coulometer.



NOTE

English is set as the default dialog language when the instrument is switched on for the first time.

Information about changing the dialog language can be found in the chapter *Selecting the dialog language* of the detailed manual.

Proceed as follows:


Press the power switch on the left-hand side of the back panel of the 917 Coulometer.

The 917 Coulometer is initialized. A system test is performed. This process takes some time.



NOTE

If a buret unit is connected, then a request appears to carry out the **Prepare** function:

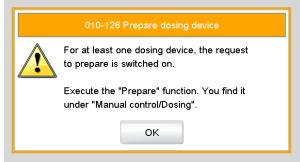


All tubings and the cylinder are rinsed with the **Prepare** function.

The preparing of the buret unit is described in the chapter *Manual control* of the detailed manual.

• Confirm the message with **[OK]**.

The main dialog is displayed:

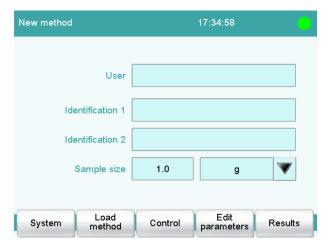

Switching off the instrument



CAUTION

The 917 Coulometer must be switched off by pressing the power switch on the rear of the instrument before the electricity supply is interrupted. If this is not done, then there is a danger of data loss.

32 ••••••

6 Operation

Proceed as follows:

Press the power switch on the left-hand side of the back panel of the 917 Coulometer.

The current data is saved and the system is shut down. This process takes just a short time. At the same time, all other instruments connected to the 917 Coulometer via a USB cable are also being switched off (except the 885 Compact Oven SC).

6.2 Fundamentals of operation

6.2.1 Touch-sensitive screen

The entire 917 Coulometer user interface is touch-sensitive. Simply touch a few of the buttons on the interface to learn how a touch-sensitive screen reacts. You can always return to the main dialog by touching [$\widehat{\mathbf{h}}$].

In order to enable an element on the 917 Coulometer user interface, just touch the screen with your fingertip, the eraser of a pencil or a stylus (special tool for operating instruments with touch-sensitive screens).



CAUTION

Never touch the touch screen with a pointed or sharp object such as a ballpoint pen.

In the default setting, the software is configured in such a way that an acoustic signal will be generated every time an active control is touched. This setting can be deactivated in the system settings.

6.2.2 Display elements and controls



The following display elements and controls are available:

Table 1 Fixed keys which are always available

	[Home] always opens the main dialog.
\(\(\)	[Back] saves the entry and opens the next-higher dialog page.
?	[Help] opens the online help for the dialog displayed.
	[Print] opens the printing dialog.
₹	[Manual] opens the manual control.
	[Stop] cancels the running determination.
\triangleright	[Start] starts a determination.

The file name of the loaded method, the time and the system status are displayed in the main dialog in the **Title bar**.

In the other dialogs, the title bar shows the headings of the next upper level and of the displayed dialog. This is an aid for orientation during navigation through the user dialog.

Table 2 Screen elements



Buttons open a new dialog when they are tapped.

34 •••••• 917 Coulometer


-----6 Operation

Delete	Inactive buttons with gray lettering indicate that the respective function is not available at the moment.
1.0	Input fields open an input dialog when tapped.
	Tapping on the selection symbol opens a selection list.
	A check box can also be activated or deactivated by tapping on it.

Status display 6.2.3

The current status of the system is displayed in the upper right-hand corner of the title bar.

Table 3

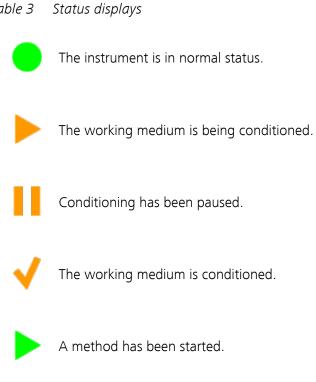



A method has been paused.

6.2.4 Entering text and numbers

In the editing dialog for text or numerical input, enter the individual characters by tapping in the input field. The following functions are available:

Text editor



Table 4 Editing functions

Editing function	Description
[OK]	The modification is applied and the editing dialog is exited.
[Cancel]	The editing dialog is exited without applying the modification.
[Delete entry]	The content of the input field is deleted completely.
[0]	The character in front of the cursor is deleted.
[⇔]	The cursor within the input field is shifted to the left by one character.
[⇔]	The cursor within the input field is shifted to the right by one character.
[az]	The lower-case letters are displayed. The label changes to [AZ] . The upper-case letters are displayed again by tapping.
[0-9]	Numbers and mathematical characters are displayed.
[Special charac- ters]	Special characters are displayed. You can use the button [More] to navigate through all available characters.

36 ••••••

6 Operation

Number editor

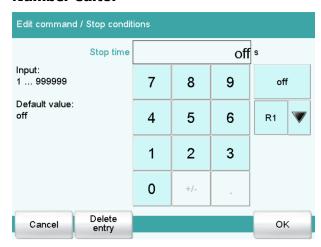


Table 5 Editing functions

Editing function	Description
[OK]	The modification is applied and the editing dialog is exited.
[Cancel]	The editing dialog is exited without applying the modification.
[Delete entry]	The content of the input field is deleted completely.
[off]	If not only numbers but also special values (e.g., off) can be entered, then the corresponding buttons will be shown to the right of the numerical keypad.
[R1]	For many parameters, a result previously defined in the method can also be entered in place of a number. You can select the result variable by touching [R1] .



NOTE

A commercially available USB keyboard can be connected to make text and numerical input easier.

7 Maintenance

Maintenance

The electronic and mechanical functional groups of Metrohm instruments can and should be checked by specialist personnel from Metrohm as part of a regular preventive maintenance schedule. Please ask your local Metrohm representative regarding the precise terms and conditions involved in concluding a corresponding maintenance agreement.

For detailed information on this topic, please visit www.metrohm.com.

Cleaning

Cleaning the surfaces of the instrument

Prerequisites

- The instrument is disconnected from the power grid.
 - 1 Clean the surfaces with a damp cloth.



NOTE

Water or ethanol can be used as a cleaning medium.



NOTE

The connectors at the rear of the instrument must only be cleaned with a dry cloth.

38 ••••••

Index

Index

800 Dosino	24	Manual 9	Power supply unit	
805 Dosimat	24	Print 9	Connect	17
		START 9	Power switch	10
Α		STOP 9	Printer	26
Addition and aspiration tube				
Mount	16	G	R	
Adsorber tube		Generator electrode 10, 18	Remote Box	
Fill	12	Fill 15	Connect	25
В		<u>I</u>	S	
Balance	27	Indicator electrode 10, 20	Safety	5
Barcode reader		Instrument	Safety instructions	6
Connect	28	Switch off 32	Sensor	
BRC	2	Switch on 31	Connect	18
			Service	6
C		K	Stirrer	
Connect		Keyboard	Connect	25
Balance	27	Connect 28	Supply voltage	7
Barcode reader	28	KFC 2	Switch off	
Dosing device	24		Switch on	31
MSB devices		M		
PC keyboard		Metrohm Serial Bus 10	Т	
Printer		Metrohm Serial Bus MSB, see also	Temperature sensor	. 10, 22
Remote Box	25	"MSB" 23	Connect	
Stirrer or titration stand		Molecular sieve	Text input	36
USB devices	26	Replace 13	Titration cell	
USB hub		MSB	Equip	14
		Connect devices 23	Fill	
D		MSB connector 10	Insert	
Display	9		Set up	12
Dosing device		N	Titration mode	
Connect	24	Network connection 30	BRC	2
		Numerical input	KFC	2
E			Titration stand	
Electrode		0	Connect	
Connect	18	Operation	Touch screen	
Electrostatic charge		General 33	Type plate	
Ethernet connector 1			.,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,	
	,	P	U	
F		PC keyboard	USB	
Fixed key		Connect 28	Connect devices	26
Back	9	Power connection 10	USB connector	
Help		Power supply	USB hub	
Home		Power supply unit 17	Connect	26
				5