

1 Purpose

This document describes the differences between the new software version **MagIC Net 4.2 Patch 1 (Build 152-1)** and its predecessor **MagIC Net 4.2 (Build 152)**.

2 Fixed bugs

Workplace program part

- Performing a determination in MagIC Net 4.2, in which a device was controlled via RS232, caused the software to freeze sporadically.


Database program part

- If a monitored result falls below or exceeds a limit value, then the determination is highlighted in red in the overview. If this monitored result was adjusted by reprocessing so that it no longer fell below or exceeded a limit value, then the determination would nevertheless remain highlighted in red if "Reprocessing with history" was deactivated ("off") in the security settings.
- When exporting with SLK file format, an error message appeared, despite the successful export of the data to the SLK file. The exported values were then displayed in black in the SLK file, even though they should have been displayed in green or red by the monitoring.

3 Compliance

The current software version does not contain any modifications that affect conformity of **MagIC Net** regarding 21 CFR Part 11 and GAMP.

Herisau, August 25, 2025



P. Hunziker
Vice President,
Head of Development



Dr. M. Saba
Head of Quality Management