

## 1 Purpose

This document describes new features and fixed bugs in the following new software version in comparison with the previous version.

- New version: 4.1.1.362
- Old version: 4.1.1.329

## 2 New features

None

## 3 Improvements

None

## 4 Fixed bugs

| Instrument            | Topic                  | Description   |
|-----------------------|------------------------|---|
| All Raman instruments | Memory leaks           | Memory leaks in the dll have been fixed. They led to the application finish after 10,000 cycles.  |
| All Raman instruments | Communication recovery | If the communication to the instrument was lost, the only way to re-establish the connection was to restart the instrument. The communication is now recovered automatically by the application itself. |
| All Raman instruments | Performance test       | PVR file compatibility between Vision and BWSpec has been improved.   |
| All Raman instruments | Performance test       | Results from performance test were calculated incorrectly.  |

## 5 Known bugs

None



## 6 Compliance

The current software version does not contain any modifications that affect conformity of Vision regarding GAMP.

Requirements of 21 CFR Part 11 can only be activated when a Vision Pharma License has been used for the installation of Vision.