

1 Purpose

This document describes the innovations in the following new firmware version:

- 5.863.0030

in comparison with the firmware version:

- 5.863.0022

2 What's new

Topic	Description
New processor	The device runs with a new processor. Devices with firmware version 5.863.0020 or older can therefore not be updated to firmware version 5.863.0030. A new device with a new processor is required for using firmware version 5.863.0030.
Methods and backups	Methods and backups which were created on program version 5.863.0030 are not reverse-compatible.
USB Plug & Play	USB flash drives and printers can be connected or removed at any time. The device does not need to be switched off before a USB flash drive or a printer is connected or removed.
Languages installed	All available languages are installed on the device. No additional language files need to be loaded. The following languages are available: <ul style="list-style-type: none">▪ German▪ English▪ French▪ Spanish▪ Portuguese▪ Italian
Program files directory	New program files must be stored in the 863 ► SwUpdates directory.

3 Fixed bugs

None.

4 Known issues

Topic	Description
Restoration after backup	In firmware version 5.863.0030, if a backup from a lower firmware version is restored, then the settings under System ► External devices and System ► Lift will not be applied.
