

Release Notes

EN

1 Purpose

This document describes the fixed bugs and problems as well as improvements in the following new software versions in comparison with the previous versions.

- New: 2.0
- Old: 1.1

2 New features

Topic	Description
Support for additional instruments	Connecting XDS MasterLab Analyzer
	Connecting XDS OptiProbe Analyzer
	Connecting XDS SmartProbe Analyzer
	Connecting XDS Rapid Solids Analyzer (with and without spot size)
	Connecting XDS RapidContent Analyzer (with and without spot size)
	Connecting XDS Multivial Analyzer
Pharma compliance	Signing of measurements according to 21 CFR part 11
	Electronic signatures (Level, Reason, Login, User rights)
	Electronic signature of configuration change before publish (on instrument configuration changes)
	Electronic records of diagnostics tests
	User Login according to 21 CFR part 11
	All measurements available in the database at all time
	View of all instrument configuration changes

Topic	Description
Installation	Installation process of Vision Air Server redesigned and simplified.
Reports	All reports are by default opened as PDF.
DS2500	Diagnostics test on start-up of the software will be skipped, if the instrument has not been turned of.
Server configuration	Server configuration can now be edited after an installation as well.
Tools section in Vision Air Routine	Usability of Tools section in Vision Air Routine has been improved.
Sample registration page	Usability of the sample registration page in Vision Air Routine has been improved.

The current software version does now contain modifications that allow conformity of Vision Air regarding 21 CFR Part 11 and GAMP.

Metrohm AG, Ionenstrasse 6, CH-9100 Herisau/Switzerland

Metrohm AG holds the SQS certificate ISO 9001 – Quality management system for the development, production and sale of instruments and accessories for analyses.

Herisau, March 1st, 2017



P. Hunziker
Vice President,
Head of Development

M. Saba
Head of Quality Management