

## 1 Purpose

This document describes the changes, improvements and fixed bugs in the following firmware version:

- 5.900.0043

in comparison with the firmware version:

- 5.900.0042

## 2 New functions

- The manual titration (MAT) can be controlled with the 6.2107.120 Manual Dosing Controller via the 6.2148.010 Remote Box MSB.

## 3 Improvements

The following improvements were made:

- The response of the scroll bar when scrolling the on-screen display has been improved.

## 4 Fixed bugs

The following bugs were fixed:

- When using a **[MEAS Conc]** command with standard addition, the measuring point interval was changed in the background at the start of the method. Because of this, the method status changed to modified.
- Printing out the calculation report on the Custom Q3X printer lead to an error message when large calculation formulas were being used.
- The validity of an expired variable was output incorrectly in the common variable report.
- The **[SCAN]** command produced an error if the timeout was reached with the timeout action set to **[Document message]**.

## 5 Conformity

[illegible]