

1 Purpose

This document describes the changes, improvements and fixed bugs in the following firmware version:

- 5.916.0045

In comparison with the firmware version:

- 5.916.0044

2 New functions

- None

3 Rectification

- Clarification in the System Assessment Report that the operator is also involved in the fulfillment of point 11.10 (g). The 21CFR Part 11 certificate has been adapted accordingly (summary).

4 Fixed bugs

The following bugs were fixed:

- When creating a report, the **MOVE** command caused unnecessary line breaks in some settings.
- After a device restart, a report loss could occur when several PDF file reports were output quickly.
- In the device manager, the preceding characters ".\" could not be used when connecting to the shared memory for the user.
- In the file manager, the reagents were not saved during the backup.
- If a message was on the screen when the user logged out automatically, the password entry could not be confirmed when logging in.
- If the network connection to the shared memory was not used for a long time, this could result in an interruption in the network connection.

5 Conformity

[illegible]

This firmware version does not contain any changes which affect the conformity of the product with regards to 21 CFR Part 11 and GAMP.