

1 Purpose

This document describes the changes, improvements and fixed bugs and problems in the following firmware version:

- 5.917.0021

in comparison with the firmware version:

- 5.917.0020

2 Improvements

The following improvements were made:

- The maximum number of users was increased from 99 to 200.
- A backup (including Audit Trail) can be saved on a PC via the network, thereby preventing unauthorized access.
From there, the Audit Trail can be viewed directly in the Audit Trail Viewer on the PC.

3 Fixed bugs and problems

The following bugs and problems were fixed:

- While printing with the thermal printer NEO's, a connection loss could occur when weighing data was transferred with the USB/RS-232 converter for 900 Touch Control (6.2148.050).
- Some of the Chinese dialog texts were wrong.
- The monitoring of the cell constant calibration validity for conductivity sensors did not work.
- Exchange units and dosing units with corroded pins were not monitored correctly, no error messages were displayed.
- When a manual titration was carried out with the MAT command, freezing of the dialog text could occur.

4 Conformity

This firmware version does not contain any changes which affect the conformity of the product with regards to 21 CFR Part 11 and GAMP.