

## 1 Purpose

This document describes the changes, improvements and fixed bugs in the following firmware version:

- 5.917.0024

in comparison with the firmware version:

- 5.917.0023

## 2 New functions

- None

## 3 Improvements

The following improvements were made:

- LAN connection: For saving on shared memory and for printing with the network printer, the SMB protocol version 2 and 3 is supported.
- The behavior of the login dialog when the password has expired and when the maximum number of login attempts has been reached was improved. The validity of the password is only checked when logging in. If the password has expired, the user can use the **Change password** command directly.

## 4 Fixed bugs

The following bugs were fixed:

- The user interface sometimes froze when printing a report on the network printer.
- If a Remote Box and a dosing device 800/805 were connected to the same MSB connector, the method run was sometimes interrupted during the titration.
- Method with **REQUEST** command with setting Stop run **Off**: If the method was started directly from the help dialog, the **REQUEST** dialog was overwritten with the live curve and the measured values in the method run.



- **KFC** command with conditioning: If the titration was started in the status **Conditioning not OK** and if the request message was not confirmed for a longer time, the user interface responded very slowly after the message had been confirmed.

## 5 Conformity

This firmware version does not contain any changes which affect the conformity of the product with regards to 21 CFR Part 11 and GAMP.