


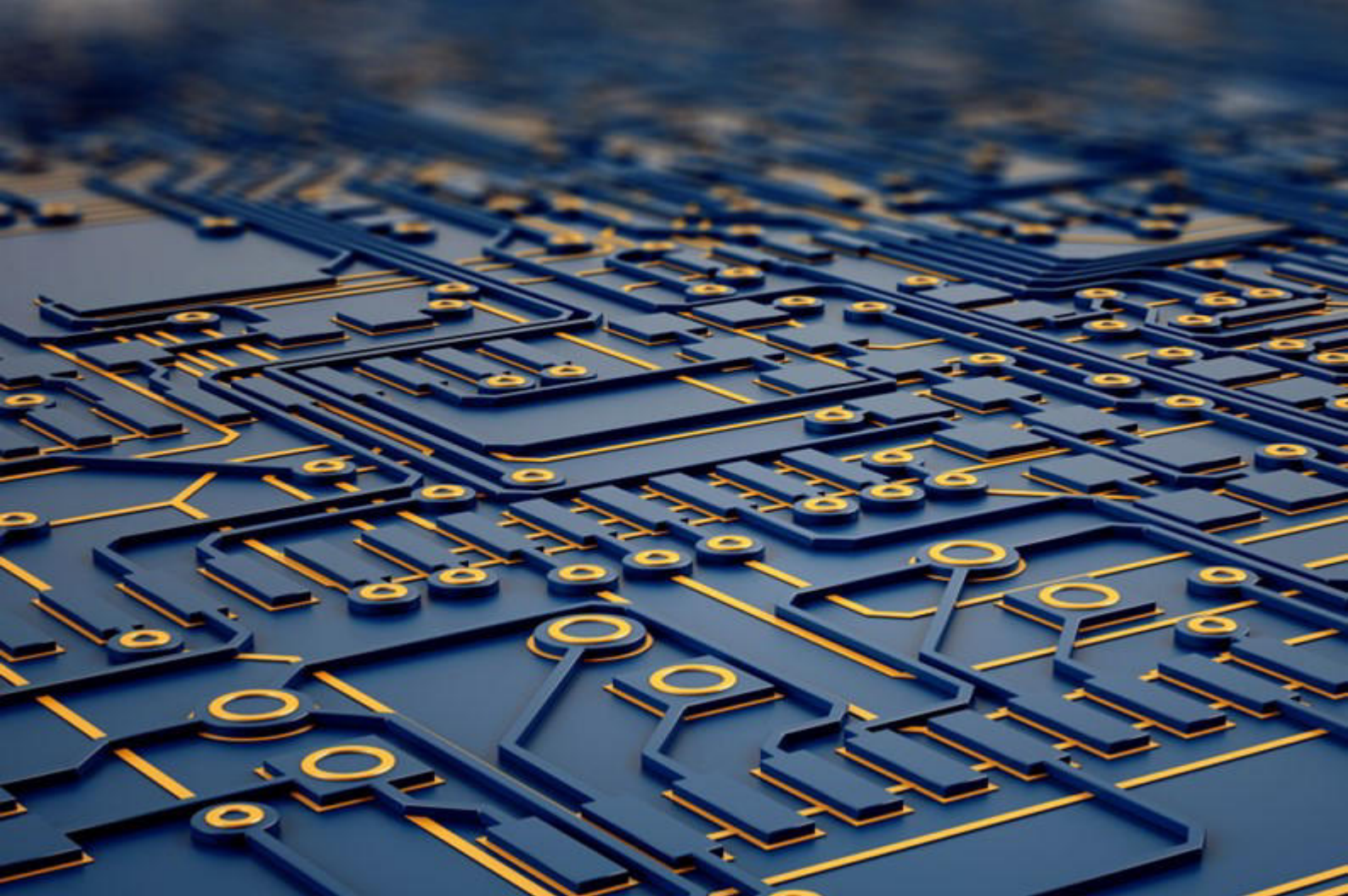

Application Note AN-V-238

無電解ニッケルめっき浴中のヒスマス (Bi) 安定剤の定量

scTRACE Gold

ENIG()ENEPIG()(Bi)
(ASV)scTRACE GoldscTRACE Gold884

Professional VA

1884 Professional VA(1)2
scTRACE Gold




1. 884 Professional VA

1.

	DP – Differential Pulse
	-0.1 V
	30 s
	0.0 V
	0.3 V
	0.15 V

- scTRACE Gold

1 mg/L30100 g/LscTRACE Gold

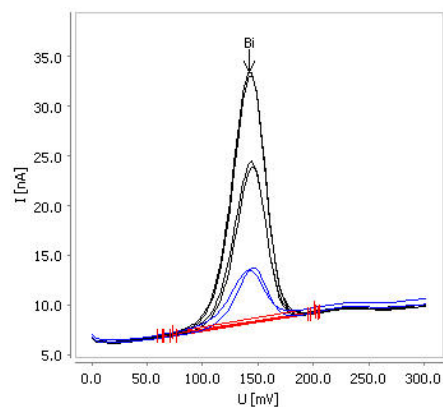

2. 1 mg/L3025 μ L10 mL

2.

Bi [mg/L]	
1 mg/L Bi	1.07

CONTACT

143-0006 6-1-1
null 9

metrohm.jp@metrohm.jp



(MME) 884 Professional VA manual
 (MME) 884 Professional VA manual proscTRACE
 Gold / viva Metrohm
 CVS ()CPVS ()CP ()
viva
 884 Professional VA manual MME pro **viva**