



Application Note AN-K-070

- ASTM D6304

Fully automated determination according to ASTM D6304

..... :
ASTM D6304 (B)

SAMPLE AND SAMPLE PREPARATION

/

EXPERIMENTAL

'''
'''
(),



Figure 1. 874 Oven Sample Processor 851 Titrando

RESULTS

„ 1.

1 ASTM D6304

(n = 4)	/ (l)	() / (l)	()/ (%)
	27.8	1.7	6.08
	44.6	0.7	1.57
	22.9	1.1	4.63
	2830.7	6.2	0.22
	18.9	2.5	13.39
	17.0	0.6	3.41

CONCLUSION

874 Oven Sample Processor 851 Titrando, ASTM

D6304 B,, 874 Oven Sample Processor ,,

CONTACT

117702
100085

marketing@metrohm.com.cn

CONFIGURATION



851 Titrando

801

(10 µg-10 mg),,,

851 Titrando

:10 µg-200 mg

OMNISiamo, GMPGLP FDA , 21 CFR Part 11



874 Oven Sample Processor

874 Oven Sample Processor · KF

OMNIS
A WHOLE NEW LEVEL OF PERFORMANCE

OMNIS 2

OMNIS Software Windows™

:

- OMNIS

-

-